


DataMate Guidebook

Data Collection using LabPro[®], Vernier Sensors, and Texas Instruments graphing calculators

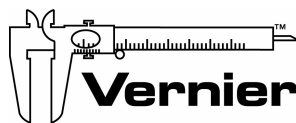

Vernier Software & Technology

13979 S.W. Millikan Way • Beaverton, OR 97005-2886

Toll Free (888) 837-6437 • (503) 277-2299 • FAX (503) 277-2440

info@vernier.com • www.vernier.com

DataMate Guidebook © 2000 Vernier Software & Technology. All rights reserved. You are entitled to reproduce parts of this book for use in your school only. No part of this book may be used or reproduced in any other manner without written permission of Vernier Software except in the case of brief quotations embodied in critical articles or reviews.

The terms Vernier LabPro and Logger *Pro* are registered trademarks of Vernier Software & Technology in the United States of America. Graphical Analysis is copyrighted by Vernier Software & Technology. CBL, CBL 2, TI Connect, and TI-GRAPH LINK are trademarks of Texas Instruments Incorporated. The term Macintosh is a registered trademark of Apple Computer, Inc. Windows is registered trademark of Microsoft Corporation.

Published by
Vernier Software & Technology
13979 SW Millikan Way
Beaverton, Oregon 97005-2886
(503) 277-2299
FAX (503) 277-2440
www.vernier.com
info@vernier.com

Third Edition 2004
First Printing
Printed in the United States of America

DataMate Guidebook

Table of Contents

DataMate Guidebook	3
Introduction	3
Required Equipment and Supported Calculators	3
Connect the Equipment and Load DataMate onto your Calculator	4
Connect the Hardware	4
Provide Power to LabPro.....	4
reset calculator (optional but recommended).....	4
Calculator Program Installation	5
TI-82, and TI-83.....	5
TI-73, TI-73 Explorer, TI-83 Plus, TI-83 Plus SE, TI-84 Plus, TI-84 Plus Silver Edition	5
TI-86	5
TI-89, TI-89 Titanium, TI-92, TI-92 Plus, and Voyage 200.....	5
LabPro Tutorials	6
Overview	6
LabPro Tutorial #1: Temperature Measurement with DataMate and Auto-ID Sensors.....	7
TI-82, and TI-83 Calculators.....	7
TI-73, TI-73 Explorer, TI-83 Plus, TI-83 Plus Silver Edition, TI-84 Plus, TI-84 Plus Silver Edition Calculators.....	7
TI-86 Calculators	7
TI-89, TI-89 Titanium, TI-92, TI-92 Plus, and Voyage 200.....	7
LabPro Tutorial #2: Using a Non Auto-ID Sensor	9
LabPro Tutorial #3: Events with Entry and Boyle’s Law	11
LabPro Tutorial #4: Collect Motion Data.....	13
LabPro Tutorial #5: LabPro as a Stand Alone Device.....	15
The DataMate Program in Detail.....	17
Overview	17
Quick Start.....	17
TI-82, and TI-83 Calculators.....	17
TI-73, TI-73 Explorer TI-83 Plus, TI-83 Plus Silver Edition, TI-84 Plus, TI-84 Plus Silver Edition Calculators.....	17
TI-86 Calculators	17
TI-89, TI-89 Titanium, TI-92, TI-92 Plus, and Voyage 200.....	17
Setup Option.....	19
Sensors	19
Data Collection Modes.....	20
Calibration.....	22
Zero	23
Save/Load.....	23
Start Option	24
Graph Option.....	25
Analyze Option.....	26
Tools Option.....	27
Quit Option.....	28
Supplemental Programs.....	29
Load DataDir to Your Calculator	29

Start the DataDir Program	30
Store a Program.....	31
Retrieve a Program from Storage	31
Delete a Program from Storage	32
Check Memory	32
Collect Garbage.....	33
Exit the Program.....	33
Supported Program Types	34
Using LabPro with Programs for the Original CBL.....	35
Sending Graphs and Data to Your Computer.....	36
Using Logger Pro and Graphical Analysis Software	36
Importing procedures for TI-73, TI-73 Explorer, TI-83, TI-83 Plus, TI-83 Plus SE, TI-84 Plus, TI-84 Plus SE, TI-86, TI-89, TI-89 Titanium, TI-92, TI-92 Plus, Voyage 200:.....	37
Appendix A: Battery and Adapter Information	39
Operating Power Requirements	39
When to Replace Batteries	39
Recommended Batteries	39
Installing the AA Batteries	39
Connecting an Optional AC Adapter.....	39
Approved AC Power Adapters	40
Building an External Battery Adapter Cable	40
Connecting an External 6-Volt Battery.....	40
Appendix B: DataMate and Calculator Specifics	41
Calculator Differences	41
Appendix C: LabPro Products, Calculators, and Sensors	42
Sensors for the LabPro and CBL 2	42
Do I Need To Purchase an Adapter?	43
Appendix D: Support and Service Information	44
International Orders.....	44
Product Service.....	44
Other Services	44
Warranty Information	44
Appendix E: DataMate Troubleshooting Guide.....	45
Appendix F: DataMate Sensor Setup Default Settings	49

Introduction

This guidebook provides step-by-step instructions for collecting real world data using LabPro¹—a portable, handheld data collection device—with a Texas Instruments graphing calculator and Vernier sensors². DataMate supports all analog sensors and the Motion Detector. (Most users never need another program. For special probes, such as photogates, rotary motion sensors, radiation monitors, and drop counters, or special features, such as graph matching, we also have other calculator programs. Installing and using supplemental programs is explained in the section titled Supplemental Programs.)

By following this guidebook, you will:

- identify the necessary equipment
- connect the equipment
- load the data collection program, DataMate, on your calculator
- collect data by following the tutorials

REQUIRED EQUIPMENT AND SUPPORTED CALCULATORS

To collect data with LabPro and a calculator, the following materials are needed:

- LabPro System: LabPro interface, AC power supply, calculator link cable, DataMate calculator program, calculator cradle, user's manual, and voltage sensor.
- A TI-73, TI-73 Explorer, TI-82, TI-83, TI-83 Plus, TI-83 Plus Silver Edition, TI-84 Plus, TI-84 Plus Silver Edition, TI-86, TI-89, TI-89 Titanium, TI-92, TI-92 Plus, or Voyage 200 graphing calculator.

Because of their Flash ROM capabilities, we recommend the TI-73, TI-73 Explorer, TI-83 Plus, TI-83 Plus Silver Edition, TI-84 Plus, TI-84 Plus Silver Edition, TI-89, TI-89 Titanium, or Voyage 200 calculators. If you are using version 2 of the AMS software for the TI-89 or TI-92 Plus, you must update to version 2.04 or newer.

- Sensors: Sensors can be divided into two basic types—analogue and digital. Examples of analogue sensors are temperature probes, pH sensors, force sensors, etc. Up to four analogue sensors can be connected to LabPro. An example of a digital sensor is a motion detector. Up to two digital sensors can be connected to LabPro.



Optional (but recommended)

- Four AA batteries for remote data collection.
- Graphical Analysis, Logger Pro, TI Connect™, and/or TI-GRAPH LINK™ software.


¹ LabPro can also be used as a computer interface using Logger Pro® software. Further LabPro resources, the LabPro Manual and LabPro Technical Reference Guide, are available on the LabPro CD.

² The CBL 2™ can also be used with Vernier sensors to collect data with Texas Instruments graphing calculators. The directions contained herein apply to both interfaces, except where noted.

Connect the Equipment and Load DataMate onto your Calculator

CONNECT THE HARDWARE

1. Insert the upper end of the calculator into the cradle.
2. Press down on the lower end of the calculator until it snaps into place.
3. Slide the bottom of the cradle onto the top of LabPro until it clicks in place.
4. Plug one end of the LabPro-to-calculator cable into the I/O port in the lower end of the calculator, and the other end into LabPro.

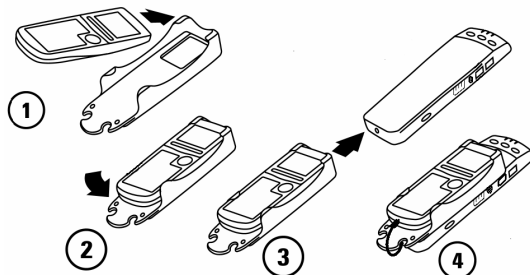

PROVIDE POWER TO LABPRO

Batteries or AC power can power LabPro. To use LabPro with batteries, remove the battery compartment cover from the underside of the interface. Insert four alkaline AA batteries and replace the compartment cover. Shortly after connecting to power, the interface will run through a self-test. You will notice a series of beeps and blinking lights (red, yellow, then green) indicating a successful power up. If the self-test is not successful, make sure that the batteries are installed correctly and that they are still useable.

To use LabPro with the AC power supply, plug the round plug on the 6-volt power supply into the side of the interface. Shortly after plugging the power supply into the outlet, the interface will run through a self-test. You will notice a series of beeps and blinking lights (red, yellow, then green) indicating a successful power up.

RESET CALCULATOR (OPTIONAL BUT RECOMMENDED)

The TI-82, TI-83, and TI-92 calculators are very limited in memory. To allow room for data storage and processing, you *must* remove all other programs and data from the calculator. Other calculators have sufficient memory to hold other data and programs along with DataMate, but to maximize the amount of data you can collect you should remove other programs. The easiest way to do this is to reset the calculator. **Warning: Resetting the calculator will remove all programs and data. Back up any needed programs or data before proceeding.**

TI-73, 73 Explorer: Press MEM. Select Reset, and then select All RAM. Select Reset.

TI-82: Press MEM. Select Reset, and then Reset again. You may need to increase screen contrast by briefly pressing and then holding down the key until the screen is legible.

TI-83: Press MEM. Select Reset, and then select All Memory. Select Reset. You may need to increase screen contrast by briefly pressing and then holding down the key until the screen is legible.

TI-83 Plus, 83 Plus Silver Edition, 84 Plus, 84 Plus Silver Edition: Press $\boxed{2\text{nd}}$ MEM. Select Reset, and then press $\boxed{\blacktriangleright}$ twice. Select All Memory. Select Reset.

TI-86: Press $\boxed{2\text{nd}}$ MEM. Select RESET. Select ALL. Select YES.

TI-89, 89 Titanium, 92, 92 Plus, and Voyage 200: Press $\boxed{2\text{nd}}$ MEM. Press $\boxed{F1}$. Select All Memory (All on the TI-92), and press $\boxed{\text{ENTER}}$. You may need to adjust screen contrast by holding down $\boxed{\blacklozenge}$ and then pressing the $\boxed{-}$ key until the screen is legible.

CALCULATOR PROGRAM INSTALLATION

When it is time to collect data, you will run a program called DataMate on the calculator. New calculators do not contain the program. You must install it, but the process is very simple. The LabPro automatically determines which kind of calculator is connected and transmits the appropriate version of DataMate.

1. Prepare the calculator to receive the program.

TI-73, TI-73 Explorer, TI-82, TI-83, TI-83 Plus, TI-83 Plus Silver Edition, TI-84 Plus, TI-84 Plus Silver Edition

Turn on your calculator and press $\boxed{2\text{nd}}$ [LINK] (on the TI-73 and TI-73 Explorer, press $\boxed{\text{APPS}}$, then select Link...). Press $\boxed{\blacktriangleright}$ to RECEIVE, then press $\boxed{\text{ENTER}}$ (Waiting... appears on your screen).

TI-86

Turn on your calculator and press $\boxed{2\text{nd}}$ [LINK], then press $\boxed{F2}$ (Waiting... appears on your screen).

TI-89, TI-89 Titanium, TI-92, TI-92 Plus, and Voyage 200

Turn on your calculator. Set it to the home screen by selecting the home key.

2. Press the TRANSFER button on the LabPro. "Receiving..." followed by a list of the loaded programs (or application) is displayed on the calculator.
3. Two beeps from LabPro will indicate the transfer is complete. Press $\boxed{2\text{nd}}$ [QUIT].

TI-82, and TI-83

Press $\boxed{\text{PRGM}}$ to see DataMate and its necessary subprograms.



TI-73, TI-73 Explorer, TI-83 Plus, TI-83 Plus Silver Edition, TI-84 Plus, TI-84 Plus Silver Edition

Press $\boxed{\text{APPS}}$ to see DataMate.




TI-86

Press $\boxed{\text{PRGM}}$ and then press [NAMES] to see DataMate and its necessary subprograms.



TI-89, TI-89 Titanium, TI-92, TI-92 Plus, and Voyage 200

Press $\boxed{2\text{nd}}$ VAR-LINK to see DataMate and its necessary subprograms.

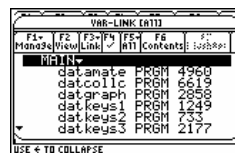

LabPro Tutorials

OVERVIEW

The tutorials will teach you how to use the DataMate program. They will explore various data collection modes using a variety of sensors. There are five tutorials that cover the most widely used forms of data collection: time graph data collection, events with entry data collection, and using the LabPro as a stand-alone device (without a calculator connected).

- DataMate automatically recognizes an auto-ID sensor, identifies the channel to which it is connected, loads a default experiment appropriate to the sensor, and shows the current reading. All active channels are shown, and the Main screen is updated as auto-ID sensors are added. In Tutorial #1 you will learn how to start the program and start collecting data right away.

Note: It is important to start with Tutorial #1, as it provides necessary setup information. Beginning with Tutorial #2 if you are using a non auto-ID sensor.

- You may have older sensors that don't have the auto-ID ability. The directions for setting up older sensors are given in Tutorial #2.
- Some experiments are not time based; instead they are event based. A pH titration is an example. Tutorial #3 will guide you through the events with entry data collection procedure using a pressure sensor.
- Tutorial #4 will show you how to set up and conduct an experiment with the Motion Detector.
- Tutorial #5 will show you how to set up LabPro as a stand-alone device for data logging, then transfer the data back to your calculator for viewing.

LABPRO TUTORIAL #1: TEMPERATURE MEASUREMENT WITH DATAMATE AND AUTO-ID SENSORS

A good way to get acquainted with LabPro is by performing a simple experiment. This particular tutorial will investigate the response time of a Temperature Probe using the DataMate program in Time Graph mode. Just follow the four easy steps and you are on your way to becoming comfortable with LabPro and DataMate.

MATERIALS

TI graphing calculator	Vernier Stainless Steel Temperature Probe
LabPro interface	One 250 mL beaker or cup
DataMate program	100 mL hot water


PROCEDURE

Connect the hardware.

1. Plug the link cable into the link port of the TI graphing calculator and LabPro.
2. Plug the Stainless-Steel Temperature Probe into Channel 1.
3. Either install four AA batteries or connect the AC adapter to LabPro.
4. Now you can start the DataMate program.

TI-82, and TI-83 Calculators

Press **PRGM** then press the calculator key for the *number* that precedes DataMate program (usually **1**). Press **ENTER**. An introductory screen will appear, followed by the main screen.

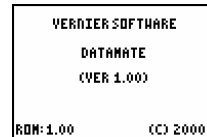

TI-73, TI-73 Explorer, TI-83 Plus, TI-83 Plus Silver Edition, TI-84 Plus, TI-84 Plus Silver Edition Calculators

Press **APPS**, then press the calculator key for the *number* that precedes DataMate program. An introductory screen will appear, followed by the Main screen.

TI-86 Calculators

Press **PRGM**, press **F1** to select < NAMES >, and press a menu key to select < DATAM > (usually **F1**). You may need to press **MORE** to see < DataMate >. Press **ENTER**. An introductory screen will appear, followed by the Main screen.

TI-89, TI-89 Titanium, TI-92, TI-92 Plus, and Voyage 200

Press **2nd** [VAR-LINK]. Use the cursor pad to scroll down to “datamate”, then press **ENTER**. Press **)** to complete the open parenthesis on the entry line and press **ENTER**. An introductory screen will appear, followed by the Main screen.

5. If the calculator cannot establish a link with LabPro, this menu appears. If you choose INTERFACE, the calculator will continue to attempt linking with LabPro. If the link still cannot be established, the error message shown below will appear.



6. Check that LabPro has power, the link cable is properly connected, and that the LabPro unit is receiving power. This screen will appear until the link problem is fixed.

```

***LINK ERROR***
MAKE SURE THE
LINK CONNECTORS
ARE FIRMLY
PUSHED IN.

[ENTER]

```

7. Press **CLEAR** to reset the program. This is the Main screen that appears after the calculator and interface are communicating. In this example, DataMate automatically identified the Temperature Probe, displayed the channel it was connected to, and displays the current live reading. DataMate also sets up a default time graph experiment for auto-ID sensors. The default settings for the temperature sensor are for a 180 second time graph.


```

CH 1:TEMP(C) 22.3

MODE:TIME GRAPH-180
-----
1:SETUP      4:ANALYZE
2:START     5:TOOLS
3:GRAPH     6:QUIT


```

8. You're ready to collect data. Select START by pressing the **2** key. (During collection, the graph is scaled to the default settings). Wait about five seconds and place the Stainless Steel Temperature Probe into the cup of hot water.

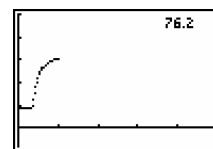

Data collection will stop after 180 seconds (or you can press the **STO▶** key to stop the data collection *before* the 180 seconds has elapsed).

9. After the data collection is complete, an autoscaled graph of the data will appear. Notice that the probe did not read the new temperature instantly; it took a little while to respond. This is a characteristic of any sensor. Use the **▶** or **◀** keys (or the cursor pad on the TI-92) to examine the data points along the displayed curve of temperature vs. time. As you move the cursor right or left, the time (X) and temperature (Y) values of each data point are displayed below the graph. Move the cursor to the point when the probe was first placed in the hot water. Record the time value. Move the cursor to find the highest temperature and record the time value. How long did it take for the Temperature Probe to stabilize at the water's temperature?

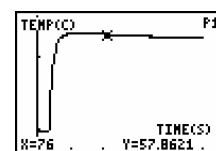

10. Press **ENTER** to return to the Main menu, then select QUIT to quit the program.

To summarize: All you need to do to collect data as a function of time with auto-ID sensors is connect the hardware, start DataMate, and select START. Time graph parameters can be changed within DataMate. Other data collection methods, using different modes, will be discussed in the tutorials that follow.

LABPRO TUTORIAL #2: USING A NON AUTO-ID SENSOR


LabPro will automatically identify auto-ID sensors; however, you may have a number of older sensors that are not automatically recognized. These will frequently be sensors with a round 5-pin DIN plug, which require a DIN-BTA adapter to connect to LabPro. The DIN-BTA adapter is a small plug with a round 5-pin DIN plug on one side, and a white, flat British Telecom plug on the other. Using these sensors is not a problem, since only a couple extra steps are required to begin data collection. This tutorial will show you how to set up non auto-ID sensors (pH is just an example) and take some data using the DataMate program in meter mode.

MATERIALS

TI graphing calculator
LabPro interface
DataMate program

distilled water

pH System with a DIN-BTA adapter.
a number of liquids in which to measure
solution (cola or any soft drink, milk,
vinegar, window cleaner, *etc.*)
beaker to hold liquids

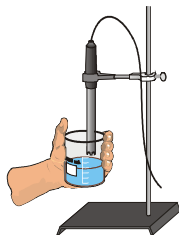

PROCEDURE

1. Connect the pH System to the DIN-BTA adapter then plug the adapter into Channel 1. Use the link cable to connect the TI graphing calculator to the interface. Firmly press in the cable ends.

```
CH 1: PH
CH 2:
CH 3:
CH 4:
DIG1:
DIG2:
MODE: TIME GRAPH-120
1:OK      3:ZERO
2:CALIBRATE
```

2. Turn on the calculator and start the DataMate program.

You will notice that DataMate has not recognized your pH System. That's because you have one of the older, non auto-ID sensors that LabPro cannot detect. It must be manually set up. Press to reset the program.

```
MODE: TIME GRAPH-20
1:SETUP   4:ANALYZE
2:START   5:TOOLS
3:GRAPH   6:QUIT
```

3. Press to select SETUP.

4. This is the Setup screen. Here you can change the sensor, units of measurement, and data collection parameters. Notice that the cursor is on CH1, indicating that you can press to modify that channel. To add a non-auto pH system to channel 1, press .

```
CH 1:
CH 2:
CH 3:
CH 4:
DIG1:
DIG2:
MODE: TIME GRAPH-20
1:OK      3:ZERO
2:CALIBRATE
```

5. A list of probes will appear. Press the number preceding PH. You will be returned to the Setup screen. By choosing a sensor type, DataMate loads the default calibration and data collection rate for that sensor. Select OK by pressing to return to the Main screen.

```
SELECT SENSOR
1:TEMPERATURE
2:PH
3:CONDUCTIVITY
4:PRESSURE
5:PID PRESSURE
6:FORCE
7:MORE
8:RETURN TO SETUP SCREEN
```

6. DataMate acts as a meter on the Main screen, and the screen updates every three seconds. The pH is displayed in the upper right-hand corner.
7. Remove the pH Sensor from the sensor storage solution bottle by unscrewing the lid. Carefully remove the bottle, leaving the cap on the sensor body. Rinse the tip of the sensor with distilled water and place the sensor tip into a beaker containing sensor soaking solution.
8. Take the pH measurement of your different liquids. Watch the readings in the corner change. Between measurements, rinse the sensor with distilled water. Switch the liquids and write down the readings. Note that DataMate does not record data in Monitor mode. Although this tutorial does not have you use the function, by selecting START you could collect pH vs. time data.

CH 1:PH	4.47
MODE: TIME GRAPH-1.00	
1:SETUP	4:ANALYZE
2:START	5:TOOLS
3:GRAPH	6:QUIT

When you are done, rinse the sensor with distilled water and return it to the sensor soaking solution. Select QUIT and exit the DataMate program.

LABPRO TUTORIAL #3: EVENTS WITH ENTRY AND BOYLE'S LAW

The Events with Entry mode collects one data point each time you press **ENTER** and then asks you to enter a corresponding value for that point. This tutorial will guide you through the setup procedure of an experiment using Events with Entry to investigate the relationship between the pressure and volume of a gas.

MATERIALS

TI graphing calculator
LabPro interface

Vernier Gas Pressure Sensor
DataMate program

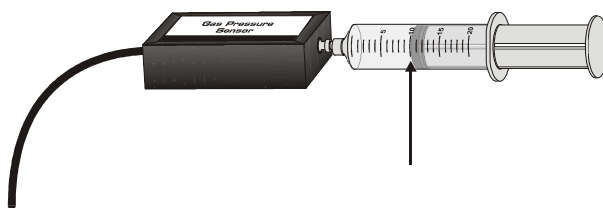


Figure 1

PROCEDURE

1. Prepare the Gas Pressure Sensor and an air sample for data collection.
 - a. Plug the Gas Pressure Sensor into Channel 1 of the LabPro. Use the link cable to connect the TI graphing calculator to the interface. Firmly press in the cable ends.
 - b. With the 20 mL syringe disconnected from the Gas Pressure Sensor, move the piston of the syringe until the front edge of the inside black ring (indicated by the arrow in Figure 1) is positioned at the 10.0 mL mark.
 - c. Attach the 20 mL syringe to the valve of the Gas Pressure Sensor. The Vernier Gas Pressure Sensors have a white stem protruding from the end of the sensor box—attach the syringe directly to the white stem with a gentle half-turn.
2. Turn on the calculator and start the DataMate program.

Notice how the program automatically identified the Pressure Sensor (an auto-ID sensor). It displays the channel it is connected to and the current reading. Press **CLEAR** to reset the program.

```
CH 1: PRESS(KPA) 102.2
MODE: TIME GRAPH-100
1: SETUP      4: ANALYZE
2: START     5: TOOLS
3: GRAPH     6: QUIT
```


Note: If you are not using an auto-ID sensor, you will have to set it up manually. Refer to Tutorial #2 for directions.

3. To collect pressure vs. volume data for Boyle's law, you need to collect the data in Events with Entry mode. The current mode is time graph, so select SETUP.
4. This is the Setup screen. Here you can change the data collection mode.

To select MODE, press **▲** once and press **ENTER**.

```
CH 1: GAS PRESSURE(KPA)
CH 2:
CH 3:
CH 4:
DIG1:
DIG2:
MODE: TIME GRAPH-100
1: OK      3: ZERO
2: CALIBRATE 4: SAVE/LOAD
```

- A list of modes will appear in the SELECT MODE screen. Press the number preceding EVENTS WITH ENTRY. You will be returned to the Setup screen. Then select OK to return to the setup screen.

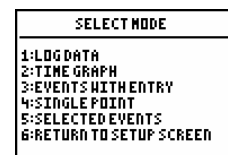

- Select OK to return to the Main screen.


You are now ready to collect pressure and volume data. It is best for one person to take care of the gas syringe and for another to operate the calculator.

- Select START. Notice that a graph is not initially displayed; instead, it is going to wait for you to collect at least two data points.

- Move the piston so the front edge of the inside black ring (indicated by the arrow) is positioned at the 5.0-mL line on the syringe.



Hold the piston firmly in this position until the pressure value displayed on the calculator screen stabilizes (in this example, the number to the right, 194.32 kPa). The number to the left indicates the sample number.

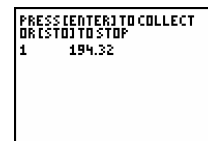

- Press **ENTER**, which tells DataMate to record the pressure. Type in “5”, the gas volume (in mL) on the calculator. Press **ENTER** to store this pressure-volume data pair.



To collect another data pair, move the syringe to 7.5 mL. When the pressure reading stabilizes, press **ENTER** and enter “7.5” as the volume.


Repeat the above procedure for volumes of 10.0, 12.5, 15.0, 17.5, and 20.0 mL.

As you work your way through the experiment, you get to view the data being collected (see the picture on the right). Notice the sample number (5), the value given (68.9 kPa) and the graph in progress.



- When you are done, press **STO** to end data collection.

- Move the cursor right or left (using the arrow keys) to examine the data pairs on the displayed graph.

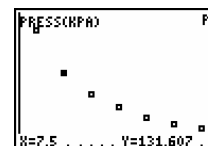

- Press **ENTER** to return to the Main menu, then select QUIT to quit the program.


LABPRO TUTORIAL #4: COLLECT MOTION DATA

The Vernier Motion Detector measures the distance to an object using echoing ultrasonic sound waves. LabPro then calculates the velocity and acceleration. This tutorial will show you how to change the length of an experiment and make a graph of a person walking.

MATERIALS

TI graphing calculator
LabPro interface

DataMate program
Motion Detector

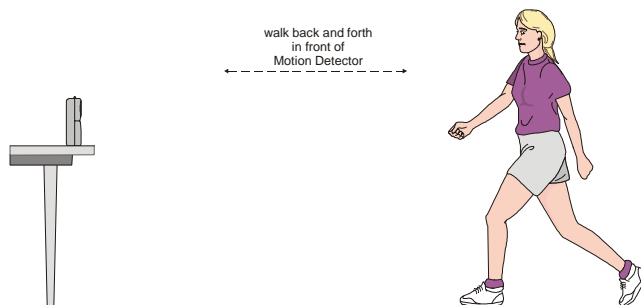

PROCEDURE


1. Place the Motion Detector so that it points toward an open space at least 4 m long.
2. Connect the Motion Detector to the DIG/SONIC 1 port of the LabPro interface. Use the black link cable to connect the interface to the calculator. Firmly press in the cable ends.
3. Turn on the calculator and start the DataMate program. Press **CLEAR** to reset the program.
4. Slowly move your hand towards and away from the Motion Detector and watch the right-hand corner of the screen as DataMate gives a live readout of the distance (in meters). Keep your hand at least 0.4 m from the detector; if you get closer than that the readings will be unreliable.
5. The system is set to collect data for five seconds, but we want a different experiment length.
 - a. Select **SETUP**.
 - b. Using the arrow keys, position the arrow next to **MODE** and press **ENTER**.
 - c. Select **TIME GRAPH**.
 - d. You will be presented with the current time settings. Select **CHANGE TIME SETTINGS**.
 - e. Type in .15 for the time interval and press **ENTER**.
 - f. Type in 100 for number of samples and press **ENTER**.
 - g. You will be presented with the new time parameters. The experiment length is now 15 seconds. Select **OK** to return to the Main screen.

```
DIG1: MOTIONCM) 1.479
MODE: TIME GRAPH-5
1: SETUP      4: ANALYZE
2: START      5: TOOLS
3: GRAPH      6: QUIT
```


```
CH 1:
CH 2:
CH 3:
CH 4:
DIG1: MOTIONCM)
DIG2:
MODE: TIME GRAPH-15
1: OK         3: ZERO
2: CALIBRATE
```

```
TIME GRAPH SETTINGS
TIME INTERVAL: .15
NUMBER OF SAMPLES: 100
EXPERIMENT LENGTH: 15
1: OK         3: ADVANCED
2: CHANGE TIME SETTINGS
```


6. Select START to collect data. You will notice a fast clicking sound coming from the Motion Detector. Wait one second, then walk slowly away from the Motion Detector. Stop for a moment, then walk toward the Motion Detector. The screen will give a live graph.

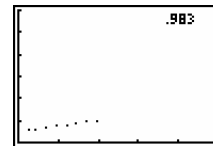

7. After data collection is complete, you will be presented with a screen asking you which graph to view. For starters, position the arrow next to DIG-DISTANCE, and press .

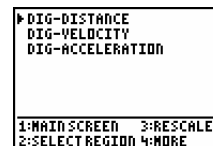


A graph (which may look something like this) of distance vs. time will appear. Press to return to the graph selection screen.

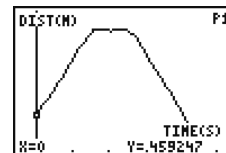

8. Press to move the cursor next to DIG-VELOCITY and press again.

A graph of velocity vs. time will appear.

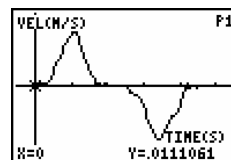

9. Press then select OK to return to the Main menu, and then select QUIT to quit the program.

LABPRO TUTORIAL #5: LABPRO AS A STAND ALONE DEVICE

The Quick Set-Up feature of LabPro lets you collect data without a calculator connected to the interface. Only auto-ID sensors can be used with Quick Set-Up.

In this mode, LabPro samples at default rates pre-set in DataMate. Data will be taken continuously and stored in the memory of LabPro. If you want to stop data collection earlier than the pre-set time, press the START/STOP button. In this experiment, you will be measuring hand temperatures of two people. LabPro will collect 99 points of data at a rate of one point every second for 99 seconds. Data collection time may be different for other sensors. Refer to Appendix F for sensor sample intervals.

EQUIPMENT

LabPro interface with batteries
DataMate program
Vernier Stainless Steel Temperature Probe

PROCEDURE


1. Disconnect LabPro from the calculator.
2. Connect the Stainless Steel Temperature Probe (an auto-ID probe) to LabPro.
3. Press [QUICK SETUP] on the LabPro. You will hear two beeps. LabPro deletes any data in memory and checks the attached probes. When the yellow light flashes, it is ready to begin collecting data.
4. Press [START/STOP]. The green light flashes to show that LabPro is collecting data.
5. Hold the tip of the probe between your fingers for thirty seconds, and then have someone else hold it.
6. When you finish collecting data, press [START/STOP] again, or wait until the 99 seconds have elapsed.
7. Transfer the data to a calculator.
 - a. Connect the calculator to LabPro with the link cable.
 - b. On the calculator, start the DataMate program or application. You will see the screen at the right.
 - c. Press [ENTER]. You will see the Main screen showing any previously setup sensors and data collection modes.
 - d. Select TOOLS.

```
DATA COLLECTION IS DONE.  
CHOOSE THE TOOLS OPTION,  
THEN CHOOSE RETRIEVE DATA.  
[ENTER]
```

```
CH 1: TEMP(C)  
  
MODE: LOG DATA  
-----  
1: SETUP      4: ANALYZE  
2: START     5: TOOLS  
3: GRAPH     6: QUIT
```

```
TOOLS  
-----  
1: STORE LATEST RUN  
2: RETRIEVE DATA  
3: CHECK BATTERY  
4: RETURN TO MAIN SCREEN
```

- e. Select RETRIEVE DATA. The data are retrieved from the memory of LabPro.

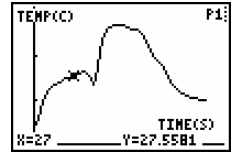

1. Scroll the cursor across the graph using the calculator's arrow keys. At which point did the Stainless Steel Temperature Probe change hands?

Note: If you start DataMate and connect a calculator during data collection, a screen showing the collection status will appear.

```
COLLECTING DATA
DIG1: 2.11243

PRESS (STOP) TO STOP.
PRESS (ENTER) TO QUIT BUT
CONTINUE COLLECTING.
```

The DataMate Program in Detail

OVERVIEW

This section describes DataMate in detail. Although the program is designed to enable you to gather data with one or two pushes of a button, there are a number of options (setting up sensors, data collection modes, etc.) that may require more detailed explanation.


We recommend that you read this section of the manual in order from start to finish.

QUICK START

The following section illustrates the fastest way to collect data using DataMate with auto-ID sensors³. Both the LabPro and CBL 2 interfaces use the DataMate program. The directions are the same for both, except as noted.

TI-82, and TI-83 Calculators

Press **PRGM** then press the calculator key for the *number* that precedes DataMate program (usually **1**). Press **ENTER**. An introductory screen will appear, followed by the Main screen.

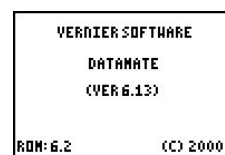

TI-73, TI-73 Explorer TI-83 Plus, TI-83 Plus Silver Edition, TI-84 Plus, TI-84 Plus Silver Edition Calculators

Press **APPS**, then press the calculator key for the *number* that precedes DataMate program. An introductory screen will appear, followed by the Main screen.


TI-86 Calculators

Press **PRGM**, press **F1** to select <NAMES>, and press a menu key to select <DATAM> (usually **F1**). You may need to press **MORE** to see <DATAM>. Press **ENTER** to go to the Main screen

TI-89, TI-89 Titanium, TI-92, TI-92 Plus, and Voyage 200

Press **2nd** [VAR-LINK]. Use the cursor pad to scroll down to “datamate”, then press **ENTER**. Press **)** to complete the open parenthesis of “datamate(” on the entry line and press **ENTER** to go to the Main screen.

If the calculator cannot establish a link with LabPro, this menu appears. If you choose INTERFACE, the calculator will continue to attempt linking with LabPro. If the link cannot be established, the error message shown below will appear.

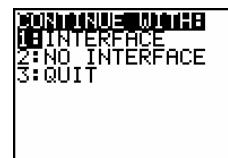

If you choose WITHOUT INTERFACE, the program will proceed to the Main screen. Under these circumstances the SETUP and START options on the Main screen will be disabled.

If you choose QUIT, the program will return you to the home screen.

³ The instructions use the DataMate program for the TI-83 and show TI-83 screen examples. DataMate programs are available for TI-73, TI-73 Explorer, TI-82, TI-83 Plus, TI-83 Plus Silver Edition, TI-84 Plus, TI-83 Plus Silver Edition, TI-86, TI-89, TI-89 Titanium, TI-92, TI-92+, and Voyage 200

If the calculator fails to communicate with LabPro, this error message appears. Check that LabPro has batteries or is powered by the AC adapter and that the link cable is properly connected. This screen will appear until the link problem is fixed.

```

***LINK ERROR***
MAKE SURE THE
LINK CONNECTORS
ARE FIRMLY
PUSHED IN.

[ENTER]
    
```

This is the Main screen. The top half of the screen shows the current sensor setup and data collection mode. The bottom half displays the menu options. A low battery icon appears in the lower right hand corner if the LabPro batteries require replacement.

```

CH 1:TEMP(C)    21.82

MODE:TIME GRAPH-100
-----
1:SETUP      4:ANALYZE
2:START     5:TOOLS
3:GRAPH     6:QUIT
    
```

The program will default to that sensor's particular Time Graph settings. In this example, you have a Temperature probe, reading in Celsius, and collecting data as a function of time for 100 seconds. The Main screen supports a meter mode displaying readings for active probes every few seconds, as displayed in the upper right-hand corner. The meter option can be toggled on and off by pressing + on the calculator.

```

CH 1:TEMP(C)    21.82

MODE:TIME GRAPH-100
-----
1:SETUP      4:ANALYZE
2:START     5:TOOLS
3:GRAPH     6:QUIT
    
```

DataMate will automatically identify auto-ID sensors, display the channel each is connected to and the current reading. Only active channels will be displayed. The screen will update as new auto-ID sensors are connected to the interface. Older sensors that do not have auto-ID need to be manually set up. Photogates, graph matching, Rotary Motion sensors, and radiation monitors require separate programs. These programs are described in the Supplemental Programs section.

You can reset the DataMate program by pressing the CLEAR button on the calculator. This reloads the calibration information onto the interface, but it also resets the DataMate program, which resets the data collection parameters. It is important to press CLEAR before setting up DataMate for a new experiment. If you do not want to lose the data collection parameters use the following method.


The second method only reloads the calibration and does not reset the DataMate program. To use this method, simply press the GRAPH button on the 73/82/83/84 or 86 calculator or the 0 button in the 89/92 or Voyage 200 while on the main screen of DataMate.

If too much time has passed while on the Main screen or on other screens, the message at right appears. If you do not respond, the calculator and LabPro will shut down to conserve batteries. If you want to continue after this happens, turn the calculator back on and select YES. When you select YES, the program resumes. If NO is selected, the program quits.

```


*****
BATTERIES
LOW
2: YES
3: NO
    
```

If you select the START option on the Main screen, LabPro will start graphing your data.

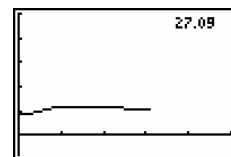

Tip: If at any time you wish to stop the data collection and view what has been collected, press the STO key.

When the timed data collection is complete, an autoscaled graph of your data is shown. Using the ← and → keys on the calculator, you can move the cursor along the graph and get readings for the x and y axes.

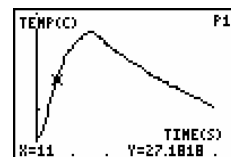

SETUP OPTION

Sensors

LabPro will automatically identify auto-ID sensors; however, you may have a number of older sensors that are not automatically recognized. These will frequently be sensors with a round 5-pin DIN plug, which requires a DIN-BTA adapter to connect to LabPro. The DIN-BTA is a small plug with a round 5-pin DIN plug on one side, and a white, flat British Telecom plug on the other. Using these sensors is not a problem, since only a couple extra steps are required to set up experiments.

NOTE: When connecting sensors to analog channels, you should use the channels in numerical order. In other words, connect the first sensor to channel 1 (CH1), the second sensor to channel 2 (CH2), and the third sensor to channel 3 (CH3). If you are using only one sensor, it should be connected to channel 1.

If the sensor is auto-ID, the channel number and sensor type automatically display on the Main screen. If necessary, change the data collection mode as described in the following section titled Data Collection Modes. Otherwise, you are ready to collect data.

If the sensor is *not* auto-ID, or if you want to change the *type* of sensor that has been auto-ID'd, follow the steps below to tell LabPro which sensor is connected.

1. On the DataMate Main screen, select SETUP.

NOTE: Here is where you will find the only difference between the LabPro and CBL 2 versions of DataMate. The LabPro setup (above) has six available channels, while CBL 2 (right) has four available channels.

- To add a non-auto-ID sensor, use the cursor keypad and position the cursor next to the channel to be set up. Press **ENTER**.
 - A list of sensors will appear.
 - Select a sensor or select MORE to see another list of sensors.
 - Selecting some sensors may produce another set of choices from which you can select a particular sensor or unit.
2. DataMate contains default time graph settings for each sensor. You can change the mode in which the collected data is graphed, and you can change the type of triggering used in the experiment. This is explained in the following section.

```
CH 1:TEMP(C)      22.3
MODE:TIME GRAPH-100
1:SETUP      4:ANALYZE
2:START      5:TOOLS
3:GRAPH      6:QUIT
```

```
▶ CH 1:STAINLESS TEMP(C)
  CH 2:
  CH 3:
  CH 4:
  DIG1:
  DIG2:
MODE:TIME GRAPH-100
1:OK          3:ZERO
2:CALIBRATE  4:SAVE/LOAD
```

```
▶ CH 1:STAINLESS TEMP(C)
  CH 2:
  CH 3:
  CH 4:
  DIG :
MODE:TIME GRAPH-100
1:OK          3:ZERO
2:CALIBRATE  4:SAVE/LOAD
```

```
▶ CH 1:STAINLESS TEMP(C)
▶ CH 2:
  CH 3:
  CH 4:
  DIG1:
  DIG2:
MODE:TIME GRAPH-100
1:OK          3:ZERO
2:CALIBRATE  4:SAVE/LOAD
```

```
SELECT SENSOR
1:TEMPERATURE
2:PH
3:CONDUCTIVITY
4:PRESSURE
5:FORCE
6:HEART RATE
7:MORE
8:RETURN TO SETUP SCREEN
```

```
TEMPERATURE
1:DIR CONNECT TEMP(C)
2:DIR CONNECT TEMP(F)
3:EXTRA LONG TEMP(C)
4:STAINLESS TEMP(C)
5:STAINLESS TEMP(F)
6:THERMOCOUPLE(C)
```

Data Collection Modes

For each Vernier sensor, DataMate loads a default experiment appropriate for the sensor. The default data collection mode for all sensors is Time Graph (collecting data points at a predetermined rate for the specific sensor). To choose another mode, or change the time graph data collection parameters, position the cursor using the arrow keys opposite MODE and press **ENTER** to see the SELECT MODE screen.

```

┌ CH 1: STAINLESS TEMP(C)
├ CH 2: LIGHT 6000(LX)
├ CH 3:
├ CH 4:
├ DIG1:
├ DIG2:
├ MODE: TIME GRAPH-20
├───┬───
├ 1:OK          3:ZERO
├ 2:CALIBRATE
└───┴───
    
```

```

┌ SELECT MODE
├ 1:LOG DATA
├ 2:TIME GRAPH
├ 3:EVENTS WITH ENTRY
├ 4:SINGLE POINT
├ 5:SELECTED EVENTS
├ 6:RETURN TO SETUP SCREEN
└───┴───
    
```

In the **LOG DATA** mode, data are collected with auto-ID sensors using Quick Setup. DataMate can then retrieve the data from the interface. Follow the directions given by this screen. See Tutorial #5 for instructions on using LabPro as a stand-alone device.

```

┌ TO COLLECT DATA, QUIT THE
├ PROGRAM, ATTACH AUTO-ID
├ SENSORS, PRESS THE QUICK
├ SETUP BUTTON, THEN PRESS
├ THE START/STOP BUTTON.
└───┴───
                                [ENTER]
    
```

The **TIME GRAPH** option is used to set the sample interval and number of data points for a timed data collection. It can also be used to set advanced features such as live graph settings and triggering.

This screen appears when you choose TIME GRAPH from the SELECT MODE screen. The default time interval between samples depends on the sensor you are using. Select CHANGE TIME SETTINGS to enter a new time interval and/or number of samples.

```

┌ TIME GRAPH SETTINGS
├ TIME INTERVAL: .14493
├ NUMBER OF SAMPLES: 138
├ EXPERIMENT LENGTH: 20
├───┬───
├ 1:OK          3:ADVANCED
├ 2:CHANGE TIME SETTINGS
└───┴───
    
```

NOTE: A time graph will be displayed live unless the sampling rate is either 0.1 seconds or smaller or larger than 270 seconds, or you have chosen to not display a live graph (see below). If you are using multiple sensors and they are identical, all of the data will be displayed live in one graph. If they are not identical, only CH 1 will be displayed live.

If you select **ADVANCED** on the TIME GRAPH SETTINGS screen, the ADVANCED TIME GRAPH SETTINGS screen appears. YMIN and YMAX refer to the “window” in which the collected data is graphed. YMIN refers to the lower boundary of the graph, and YMAX refers to the upper boundary of the graph. The YMIN and YMAX values shown on the screen are the default range of the sensor in channel 1. (This will vary depending on the sensor being used. For example, for the Stainless Steel Temperature Probe, the range is -40 to 140°C .)

```

┌ ADV. TIME GRAPH SETTINGS
├ LIVE GRAPH: TEMP(C)
├ YMIN  YMAX  VSCL
├ -20   125   25
├ TRIGGERING:NONE
├───┬───
├ 1:OK
├ 2:CHANGE GRAPH SETTINGS
├ 3:CHANGE TRIGGERING
└───┴───
    
```

Follow the steps below to change advanced time graph settings:

1. To change the window range to be graphed or turn off the live graph, select CHANGE GRAPH SETTINGS.

A list of the connected sensors appears.

```

┌ SELECT GRAPH
├ 1:CH1-TEMP(C)
├ 2:CH2-LIGHT(LX)
├ 3:NONE
└───┴───
    
```

2. Press the number beside the sensor you want, or you can choose not to display a live graph by selecting NONE.

3. To change the type of triggering, select CHANGE TRIGGERING. In the example shown, there are two types of triggering:

- LabPro will trigger the start of data collection based on a change in the data being collected. (This is called threshold triggering.)
- For MANUAL TRIGGER, the LabPro will start collecting data when the START/STOP button on LabPro is pressed.
- If you choose NONE, no special triggering will be set.

```
SELECT TRIGGERING
1:CH1-TEMP(C)
2:CH2-LIGHT(LX)
3:MANUAL TRIGGER
4:NONE
```

4. Press the number beside the type of triggering you want. If you select NONE, you will return to the ADVANCED TIME SETTINGS screen.

If you select **MANUAL TRIGGER**, the triggering option is changed and you will return to the ADVANCED TIME SETTINGS screen.

If you choose to trigger on a channel, DataMate asks you to choose the trigger type.

- INCREASING means that data collection will not start until the sensor readings (such as light intensity or temperature) increase through the selected threshold (see Step 6).
- DECREASING means that data collection will not start until sensor readings decrease through the selected threshold (see Step 6).

```
TRIGGER TYPE
1:INCREASING
2:DECREASING
```

5. Press the number beside the trigger type you want.

6. Key in the number (the threshold) at which you want data collection to begin and press **[ENTER]**. (Put in a threshold value in the units of the sensor you are using such as °C for temperature or Newtons for force.) When the values of the data being collected reach this value, LabPro will begin storing data.

```
TRIGGER
THRESHOLD: ■
```

7. Key in the number (percent) of data you want the LabPro to prestore and press **[ENTER]**. Prestore is the amount of data collected before the threshold was reached that you want to keep (10 percent, 20 percent, and so on). From the time the experiment begins until the threshold is reached, LabPro collects data in a buffer. When the threshold is reached, LabPro begins storing the data it has collected and discards the data it collected before the threshold was reached unless a prestore value is entered. The ADVANCED TIME SETTINGS screen appears.

```
PRESTORE IN
PERCENT:
```

When Manual Trigger is selected, a set of instructions will be given after you select START. After selecting START, you can leave DataMate. Then you can disconnect the calculator from the LabPro unit and use LabPro for remote data collection. Data collection will begin when the Start/Stop button on LabPro is pressed. To download data to the calculator after remote data collection, reconnect the LabPro to the calculator and restart DataMate. Follow the instructions on screen.

```
REMOTE DATA COLLECTION IS
SET UP. PRESS START/STOP
BUTTON TO COLLECT DATA.
PLEASE WAIT UNTIL
DATA COLLECTION IS DONE.
CHOOSE THE TOOLS OPTION,
THEN CHOOSE RETRIEVE DATA.
[ENTER]
```

EVENTS WITH ENTRY is used for non time-based data collection such as triggering for titrations, Boyle's law, etc. Data are collected one point at a time and only when the **ENTER** button on the calculator is pressed. You are then prompted to enter a corresponding value.

SINGLE POINT data collection is used to collect one averaged data point. The interface will collect data for 10 seconds and report a single, averaged value for each sensor on the calculator screen.

SELECTED EVENTS mode collects data from each channel in response to pressing the **ENTER** key, but the user does not enter any values. Data are plotted vs. point number.

Calibration

When a sensor is selected, DataMate automatically loads the standard calibration settings. Further calibration is usually not needed.

If you choose to calibrate a sensor, position the cursor opposite the sensor and choose **CALIBRATE** from the **SETUP** screen.

There are two ways to calibrate a sensor. The first is to perform a new calibration; the second way is to manually enter the calibration values.

Almost all Vernier sensors are calibrated with two reference points. This example describes the calibration of a pH sensor. To perform a new calibration for a pH sensor, you would need two solutions with known pH values: for example, solutions with values of 4 and 10. Here are the steps you might follow:

1. On the Setup screen, to move the cursor to the sensor you want calibrated and select **CALIBRATE**.

```

CH 1:PH
CH 2:
CH 3:
CH 4:
DIG1:
DIG2:
MODE:TIME GRAPH-180
-----
1:OK          3:ZERO
2:CALIBRATE 4:SAVE/LOAD
    
```

*Note: Not all sensors can be calibrated (Stainless Steel Temperature Probe, for example). If you select a sensor that cannot be calibrated, DataMate does not respond when you select **CALIBRATE**.*

2. The current calibration will be displayed. Select **CALIBRATE NOW**.

```


          CALIBRATION
PH
CALIBRATION:LINEAR
SLOPE      INT
-3.858     13.72
-----
1:OK
2:CALIBRATE NOW
3:MANUAL ENTRY
    
```

3. Put the pH sensor in the pH 4 solution. Wait until the voltage number on the screen stabilizes and then press **ENTER**.

```

          CALIBRATE SENSOR
MONITOR VOLTAGE, WHEN
STABLE, PRESS [ENTER].
          VALUE  VOLTAGE
POINT 1:         2.779
POINT 2:
    
```


4. Key in 4 and press **ENTER**.

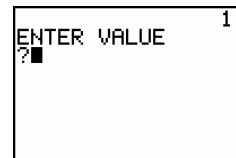

5. Repeat steps 3 and 4 for the pH-10 solution.

6. Select OK to return to the Setup screen.

To calibrate the pH sensor by entering values manually, follow these steps:

1. On the Setup screen, move the cursor to the sensor you want calibrated and select CALIBRATE.




2. Select MANUAL ENTRY.

3. Key in the slope and press **ENTER**.



4. Key in the intercept and press **ENTER**. The Calibration screen is displayed with the new values.

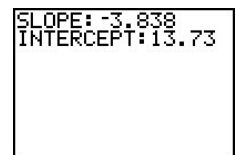

5. Select OK to return to the Setup screen.

Zero

From the Setup screen you can zero one or more sensors. After selecting the ZERO option, a screen will appear with a list of channels. Select one of the sensors you want to zero or the ALL CHANNELS option.




Another screen displaying the current reading or readings will appear. Press **ENTER** to zero the sensor or sensors.



NOTE: The ISE probes and Stainless Steel Temperature probes cannot be zeroed.

Save/Load

DataMate allows you to save experiments in LabPro's FLASH Memory, recall them later, and delete them when you no longer need them. An experiment includes all program set-ups, sensor selections, calibrations, graph settings, and so forth, as well as any data you have collected. Heart Rate Monitor and Respiration Monitor data, as well as data collected in Single Point mode, are not saved with experiments because DataMate calculates the readings.

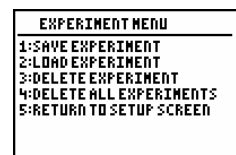

NOTE: This feature is not available for the TI-83 and TI-83 calculators.


Select **SAVE EXPERIMENT** to either save a setup with or without data. When you have entered the settings for an experiment but haven't taken data, only the settings are saved. If you have entered the settings and taken data, both are saved. Follow the steps to save an experiment.

- From the EXPERIMENT MENU screen, select SAVE EXPERIMENT.
- Type in a name and press **ENTER**. The experiment is saved, and the Setup menu reappears.



LOAD EXPERIMENT. This option allows you to call up a saved experiment to view, analyze, etc. This is useful if you want to (1) load settings so you can collect data or (2) load settings with data for analysis, etc.

- From the EXPERIMENT MENU screen, select LOAD EXPERIMENT.
- Press the number beside the experiment you want. After the experiment loads, you are taken to the main screen.

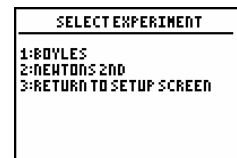


DELETE EXPERIMENT. This option deletes saved experiments one at a time.

- From the EXPERIMENT MENU screen, select DELETE EXPERIMENT.
- Press the number beside the experiment you want to delete. After the experiment is deleted, the Main menu appears.



DELETE ALL EXPERIMENTS. Not only can you delete one experiment at a time; you can also delete all the experiments that you have stored. The advantage of this feature is the time saved if several files are to be deleted. To do so, follow these steps:

- From the EXPERIMENT MENU screen, select DELETE ALL EXPERIMENTS.
- To confirm that this is what you want to do, go ahead and select YES.

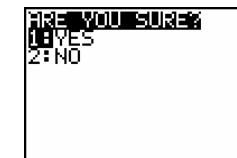

*Tip: While on the DataMate Main screen or the Setup screen, press **CLEAR** to restore DataMate's default settings, and clear previous data.*


START OPTION

Select **START** on the Main screen to begin collecting data using the existing sensor setup and data collection parameters.


GRAPH OPTION

The Graph screen automatically appears when you are finished collecting data or when Graph is chosen from the Main screen. Two conditions determine the Graph screen:

- If one analog channel is active then a graph of your data is presented when data collection is complete, or if two of the same type of analog sensors are plugged in.

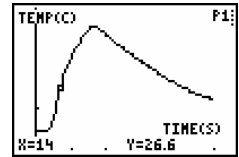

- If more than one channel is active or you have collected data using the motion detector, you will be presented with a list of available channels to view. To view a graph, select the appropriate channel from the list by using the \blacktriangle and \blacktriangledown keys on the calculator and press ENTER .

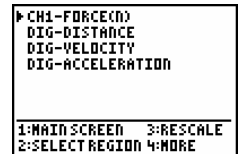


If you want to view another graph, press ENTER . The Graph screen appears again so you can choose another channel.


Selecting **MAIN SCREEN** on the Graph screen will take you back to the Main screen.

SELECT REGION on the Graph screen is used to remove unwanted data from lists, i.e. data that are outside the region of interest. This option may be necessary for curve fitting. Data outside the selected region are permanently deleted.


- Select the left bound, as viewed on the bottom of the screen, using the \blacktriangleleft and \blacktriangleright keys. When the values you want for the left bound is displayed, press ENTER .

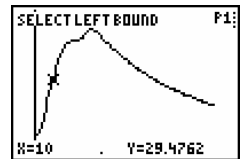

- Repeat for the right bound and press ENTER .

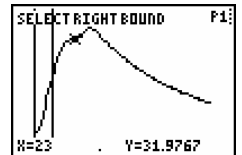


- The new graph will automatically be displayed.

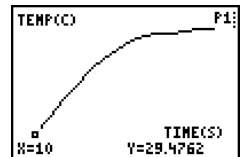

RESCALE is used to autoscale, adjust the x-axis or y-axis scaling of your graph.

- Press the number beside the scale you wish to change. You will be prompted to enter the minimum, maximum, and tic mark interval for the selected axis.

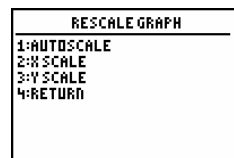

- Press **ENTER** and return to the RESCALE GRAPH screen and select another scale to see the graph with other scale options.
- When you are done viewing the graphs, press **ENTER** to return to the Rescale Graph screen, then select RETURN.

Selecting **MORE** on the Graph screen enables you to plot alternate data lists in a graph.

L1, L2, L3 and L4 refer to the lists in which your data are stored. For example, L3 vs. L1 will graph the data in List 3 versus the data in List 1.

MORE GRAPHS	
1:	L3 VS L1
2:	L3 VS L2
3:	L2 VS L3
4:	L4 VS L1
5:	L3 VS L4
6:	L2 AND L3 VS L1
7:	L2, L3 AND L4 VS L1
8:	RETURN TO GRAPH SCREEN

If multiple channels are shown, then the point, box and plus point protectors are used for lists L2, L3, and L4 respectively.

ANALYZE OPTION

The Analyze option on the Main screen produces this list of options. You can perform a curve fit, model data with a function other than curve fit functions, determine statistics over a region of data, and determine the integral over a region of data.

ANALYZE OPTIONS	
1:	RETURN TO MAIN SCREEN
2:	CURVE FIT
3:	ADD MODEL
4:	STATISTICS
5:	INTEGRAL

Selecting **CURVE FIT** from the ANALYZE OPTIONS menu produces a list of curve fit options. Selecting MORE produces another six options.

CURVE FIT	
1:	LINEAR (CH1 VS TIME)
2:	LINEAR (CH2 VS TIME)
3:	LINEAR (CH3 VS TIME)
4:	LINEAR (DIST VS TIME)
5:	LINEAR (WELD VS TIME)
6:	LINEAR (CH2 VS CH1)
7:	MORE

1:	LINEAR (CH1 VS CH2)
2:	POWER (CH1 VS TIME)
3:	POWER (CH2 VS TIME)
4:	QUAD (CH1 VS TIME)
5:	QUAD (DIST VS TIME)
6:	EXPONENT (CH1 VS TIME)
7:	PREVIOUS MENU
8:	RETURN TO ANALYZE OPT.

After a curve fit is performed, the fit parameters are displayed. Next, the fitted curve is drawn as a continuous line, and the data points are shown with box point protectors.

Note that the $t=0$ point cannot be included in power law curve fits. Use the select region option to remove the first point.

ADD MODEL graphs the collected data and overlays a graph of a general equation whose variables can be adjusted. Before choosing this option, it is necessary to quit DataMate and manually enter the general equation to be used into the Y1 equation variable. The equation may use the parameters A, B, C, D, and E as a function of X. When you select the ADD MODEL option, you will be asked to choose which graph you would like to view along with the model. The calculator will display a graph, then bring up the menu to the right.

MODEL MENU	
1:	ADJUST A
2:	ADJUST B
3:	ADJUST C
4:	ADJUST D
5:	ADJUST E
6:	RETURN TO ANALYZE MENU

The first five options will display the current value for that variable and prompt you for a new value. Each time a variable is changed, the calculator regraphs the data and the model equation. Option 6 will turn off the model and return you to the ANALYZE MENU screen. The parameters are reset upon leaving modeling.

STATISTICS finds the mean, min, max, standard deviation, and number of data points of a selected region. After selecting this option, choose a graph from the displayed list. Use the arrow keys to move the cursor and select the left bound of the region and press **[ENTER]**. A vertical line will be drawn on the graph. Now move the cursor to select the right bound and press **[ENTER]**. After the calculator has finished determining the statistics, a message will be displayed instructing you to press **[ENTER]** to continue. The numerical results will be displayed. No data are deleted in calculating statistics.

```
MEAN: 28.346
MIN: 26.000
MAX: 30.100
STD DEV: 1.322
N: 12.000

[ENTER]
```

INTEGRAL is used to integrate a section of the graph. After selecting this option, choose a graph from the list. Use the arrow keys to move the cursor and select the left bound of the region and press **[ENTER]**. A vertical line will be drawn on the graph. Use the cursor to select the right bound and press **[ENTER]**. The numerical results will be presented.

```
INTEGRAL: 312.105

[ENTER]
```

TOOLS OPTION

When Tools is selected from the Main screen, the screen to the right appears.

```
TOOLS
-----
1:STORE LATEST RUN
2:RETRIEVE DATA
3:CHECK BATTERY
4:RETURN TO MAIN SCREEN
```

After data collection, data is automatically stored in lists. For example, time is stored in List 1 (L_1) and the temperature (data) in List 2 (L_2). Subsequent data collections will overwrite those lists with the current data. You do, however, have the option to store the data (from List 2) so it won't get overwritten. Note that L_1 will not be stored. Selecting **STORE LATEST RUN** stores the previous data into List 3 so List 2 becomes free for more data. In other words, List 3 (L_3) becomes List 4 (L_4), List 2 (L_2) becomes List 3 (L_3). List 2 is then available for another run.

Note: This feature can only be used if one sensor is active in CH 1. It can only store two previous runs. The third run remains in List 2 (L_2).

RETRIEVE DATA. The Retrieve Data option is used to download data after remote data collection.

- From the TOOLS screen, select RETRIEVE DATA.
- You will get a message "Retrieving Data."
- Arrow to the data set you want to view and press **[ENTER]**. Follow the directions for viewing a graph.

```
CH1-TENPCO
CH2-PH
CH2 VS. CH1

1:MAIN SCREEN 3:RESCALE
2:SELECT REGION 4:MORE
```

CHECK BATTERY will display the status of LabPro's battery.

- From the **TOOLS** menu, press **CHECK BATTERY**.
- If battery charge is adequate you will get the following message.



QUIT OPTION

Select **Quit** to leave DataMate and return to the home screen of the calculator. On the TI-89, TI-89 Titanium, TI-92 and TI-92 Plus calculators you must press **F5** or **HOME** to return to the home screen. To see the last graph select **zoomstat** as appropriate for the calculator model.

Supplemental Programs

We have written supplemental programs for special sensors, such as photogates and radiation monitors, or for special applications, such as motion matching or managing the Flash memory of your interface.

One feature of the latest operating system for LabPro and CBL 2 is the ability to transfer the DataDir program instead of the DataMate program to your calculator from the interface. With DataDir on your calculator, you can then retrieve supplemental programs from the interface. If this is the approach that you want to take, update your interface with the most current operating system and calculator programs (available at <http://www.vernier.com/calc/flash.html>). Next reset the memory on your calculator, link the calculator and interface, and then press and hold the Transfer button on the interface. The DataDir program will be transferred to the calculator. Run the DataDir program to retrieve the supplemental programs from the interface.

Name	Use
DataGate	This program is used when you want to collect data with Photogates or Smart Pulleys™. This program will work with CBL 2 and LabPro.
DatMatch	The DatMatch program is used with the Motion Detector. The program presents distance vs. time and velocity vs. time graphs, and challenges the students to match the graph. This program will work with CBL 2 and LabPro.
DaRotary	This program is used when you want to collect data with the Vernier, Pasco or Team Labs rotary motion sensors. It supports a variety of collection modes including simultaneous analog and rotary sensor data collection. There are also a number of analysis features. This program will work with CBL 2 and LabPro.
DataRad	This program is used with the Vernier Student Radiation Monitor and our Radiation Monitor to monitor nuclear radiation. A variety of data collection modes are supported. This program will work with CBL 2 and LabPro.
FunctGen	Use a LabPro like an analog function generator.
DataDrop	Data collection with the Vernier Drop Counter.
SPECTRO	Collect data with a spectrophotometer.

The DataDir program allows you to store calculator programs in the *FLASH* memory of your LabPro/CBL 2 interface and then retrieve them later. (This is like having an “external hard drive” for your calculator.) Each LabPro and CBL 2 has approximately 600K of *FLASH* ROM available for storing experiment files and programs.

LOAD DATADIR TO YOUR CALCULATOR

There are two methods of sending the DataDir program to your Texas Instruments graphing calculator.

A. Send DataDir from a LabPro/CBL 2 Interface.

This method will only be possible if your LabPro/CBL 2 interface has operating system version 1.2 or higher. To determine the version of your operating system, connect a calculator to your LabPro/CBL 2 and start the DataMate program or one of our supplemental programs. As it starts up watch for the ROM number in the lower left of the screen. The number that appears is the version of the interface operating system.

Updates for your interface can be found on the Vernier web site at www.vernier.com/downloads/index.html and on the TI web site at education.ti.com.

1. Connect the calculator to the LabPro/CBL 2 using the short, black link cable.
2. Prepare the calculator to receive DataDir.

TI-73, TI-73 Explorer, TI-82, TI-83, TI-83 Plus TI-83 Plus Silver Edition, TI-84 Plus and TI-84 Plus Silver Edition

Turn on your calculator and press $\boxed{2nd}$ [LINK] (on the TI-73, press \boxed{APPS} , then select 1:Link). Press $\boxed{\blacktriangleright}$ to select RECEIVE, then press \boxed{ENTER} .

TI-86

Turn on your calculator and press $\boxed{2nd}$ [LINK]. Select <RECV>.

TI-89, TI-89 Titanium, TI-92, TI-92 Plus, and Voyage 200

Turn on your calculator and go to the home screen.

3. Press and hold the TRANSFER button on LabPro/CBL 2 for about 5 seconds.
“Receiving...”, followed by “DataDir”, is displayed on the calculator.
The “Done” message on the calculator and two beeps from LabPro/CBL 2 indicates the transfer is complete. Press $\boxed{2nd}$ [QUIT].

B. Send DataDir from your computer using a TI-GRAPH LINK cable.

The DataDir program is available on the TI Resource CD, on the TI web site at education.ti.com, or on the Vernier web site at www.vernier.com and can be loaded to your calculator via a TI-GRAPH LINK cable and TI-GRAPH LINK software or TI-Connect software⁵.

START THE DATADIR PROGRAM

To use the DataDir program to store and retrieve programs, your TI graphing calculator must be connected to your LabPro/CBL 2 interface.

TI-73, TI-73 Explorer, TI-82, TI-83, TI-83 Plus, TI-83 Plus Silver Edition, TI-84 Plus, TI-84 Plus Silver Edition Calculators

Press \boxed{PRGM} , then press the calculator key for the *number* that precedes DATADIR (usually $\boxed{1}$). Press \boxed{ENTER} and wait for the Main screen to load.

TI-86 Calculators

Press \boxed{PRGM} , press $\boxed{F1}$ to select <NAMES>, and press the menu key that represents DataDir. (<DATAD> is usually $\boxed{F1}$.) Press \boxed{ENTER} , and wait for the Main screen to load.

TI-89, TI-89 Titanium, TI-92, TI-92 Plus, and Voyage 200

Press $\boxed{2nd}$ [VAR-LINK]. Use the cursor pad to scroll down to “datadir”, then press \boxed{ENTER} . Press $\boxed{)}$ to complete the open parenthesis that follows “datadir” on the entry line and press \boxed{ENTER} . Wait for the Main screen to load.

An introductory screen appears briefly, then the main menu screen displays.

MAIN MENU	
1:	LOAD A PROGRAM
2:	STORE PROGRAMS
3:	DELETE A PROGRAM
4:	DELETE ALL PROGRAMS
5:	CHECK MEMORY
6:	COLLECT GARBAGE
7:	QUIT

⁵ For help using TI-GRAPH LINK software, consult *CBL Made Easy*, available on the Vernier web site at www.vernier.com/cbl/easy.html. For help using TI-Connect software see education.ti.com.

Store a Program

The program(s) you want to store on your LabPro/CBL 2 must be on your calculator, either by sending them from another calculator or transferring them from a computer using the TI-GRAPH LINK cable and the TI-GRAPH LINK software or TI-Connect software. With DataDir you can store one or more programs from your calculator to your LabPro/CBL 2 at the same time. Follow the instructions below. The key presses and screenshots in this section refer specifically to the TI-83 Plus calculator. Although the steps will be similar if you are using a different type of calculator, be sure to read and closely follow the instructions shown on your own calculator screen.

1. On the Main screen, select STORE PROGRAM. The screen to the right will appear.

```
STORE PROGRAM(S)
PRESS 2ND LINK, THEN
CHOOSE PRGM. SELECT
THE PROGRAMS TO STORE,
THEN CHOOSE TRANSMIT.
COMPLETE THIS IN
1 MINUTE.
```

2. Press [LINK].

```
SEND RECEIVE
1: All+...
2: All-...
3: Prgm...
4: List...
5: Lists to TI82...
6: GOB...
7: Pic...
```

3. Press Prgm to select a program.

```
SELECT TRANSMIT
▶ DATA DIR PRGM
  DATSHARE PRGM
  DCUALARM PRGM
  DCUAVOID PRGM
  PHOERIVS PRGM
  START PRGM
```

4. Press to move the cursor to the program you want to store and then press . A dot appears beside the program name. Repeat this step until you have selected all the programs you want to store.

```
SELECT TRANSMIT
▶ DATA DIR PRGM
  . DATSHARE PRGM
  . DCUALARM PRGM
  . DCUAVOID PRGM
  . PHOERIVS PRGM
  . START PRGM
```

5. Press to highlight TRANSMIT and press . When the program(s) have been stored, the calculator displays the message Done.

```
DATA SHARE PRGM
DCUALARM PRGM
DCUAVOID PRGM
PHOERIVS PRGM
▶ START PRGM
  Done
```

Note: The calculator exits the DataDir program to make the transfer. The transfer needs to be completed within one minute. Run the DataDir program again to see the results of the transfer.

Retrieve a Program from Storage

The DataDir program allows you to retrieve a program from storage on LabPro/CBL 2. The following instructions will guide you through this task. The key presses and screenshots in this section refer specifically to the TI-83 Plus calculator. Although the steps will be similar if you are using a different type of calculator, be sure to read and closely follow the instructions shown on your own calculator screen.

1. On the Main screen, select LOAD A PROGRAM.

```
LOAD A PROGRAM
-----
1:*DATMATCH.BXF
2:*DATAGATE.BXF
3:*DATARAD.BXF
4:*DAROTARY.BXF
5:PHDERIVS.BXF
6:DATSHARE.BXF
7:NEXT
```

2. Press the number beside the program you want to load.

3. Press [LINK].

```
PRESS 2ND LINK,
CHOOSE RECEIVE AND
PRESS ENTER].PRESS
TRANSFER BUTTON ON
INTERFACE. COMPLETE
THIS IN 1 MINUTE.
```

4. Press to highlight RECEIVE and press .

```
RECEIVE
1: All+...
2: All-...
3: Prog...
4: List...
5: Lists to TI82...
6: GDB...
7: Pic...
```

5. Press the TRANSFER button on LabPro/CBL 2.

6. When the program has been loaded on the calculator, the calculator says DONE.

Delete a Program from Storage

The DataDir program provides two options for deleting programs from storage. You can delete a single program or you can delete all the programs stored on LabPro/CBL 2.

You can only delete those programs that you have previously stored using DataDir. Deleting all programs will NOT delete DataMate.

Follow these instructions to delete a program that you have stored on LabPro/CBL 2:

1. On the Main screen, select DELETE A PROGRAM.

```
DELETE A PROGRAM
-----
1:DATSHARE.BXF
2:DCUALARM.BXF
3:DCUAVOID.BXF
4:PHDERIVS.BXF
5:START.BXF
6:RETURN TO PREVIOUS MENU
```

2. Press the number beside the program you want to delete. Only programs that you have stored on LabPro/CBL 2 using DataDir are available to be deleted. The Main screen appears.

Follow these instructions to delete ALL the programs that you have stored on LabPro/CBL 2:

1. On the Main screen, select DELETE ALL PROGRAMS.
2. The programs that you have stored are deleted and the Main screen appears.

Check Memory

The DataDir program also allows you to check available memory on LabPro/CBL 2.

Follow the directions below to check memory:

1. On the Main screen select CHECK MEMORY.



NUMBER OF AVAILABLE BYTES
IN ARCHIVE:
588256
[ENTER]

2. When you finish viewing the screen, press . The Main Screen appears.

Collect Garbage

As with most types of digital storage devices, repeated deleting and writing can lead to inefficient use of memory space. The DataDir program has a feature allowing you to clean up memory space on LabPro/CBL 2.

Follow the directions below to collect garbage:

1. On the Directory main menu select COLLECT GARBAGE.



COLLECTING GARBAGE,
PLEASE WAIT.

2. Depending on the state of LabPro/CBL 2 memory, garbage collection can take up to 5 minutes. When garbage collection is complete, the Directory main menu appears.

Exit the Program

On the Directory main menu, select QUIT. The calculator displays the message Done.

Supported Program Types

LabPro/CBL 2 and the DataDir program can handle the following calculator variable types for the listed calculators:

TI-73, TI-73 Explorer:

Store and retrieve programs. (*.73P)

Store and retrieve applications. (*.73K)

TI-82:

Retrieve programs. (*.82P)

(Does not store programs)

TI-83:

Store and retrieve programs. (*.83P)

TI-83 Plus, TI-83 Plus Silver Edition, TI-84 Plus and TI-84 Plus Silver Edition::

Store and retrieve programs. (*.8XP)

Store and retrieve applications. (*.8XK)

TI-86:

Store and retrieve programs. (*.86P)

TI-92:

Store and retrieve programs. (*.92P)

TI-89, TI-89 Titanium, and TI-92 Plus:

Store and retrieve programs. (*.89P and *.9XP)

Store and retrieve applications. (*.89K and *.9XK)

Store and retrieve assembly programs. (*.89Z and *.9XZ)

Using LabPro with Programs for the Original CBL

LabPro works with most of the existing CBL programs. Before you have your students use an older program with LabPro, test it thoroughly. If you discover problems, contact Vernier for possible updates.

There are updated versions of PHYSICS, CHEMBIO and PHYSCI for use with LabPro. They may be downloaded from the Vernier web site <http://www.vernier.com/downloads/index.html>. Using these updated programs will allow a mixture of LabPros and CBLs to be used in a classroom with a single set of student instructions.


For more information on legacy CBL programs, we recommend that you download the manual CBL Made Easy from the Vernier web site: <http://www.vernier.com/legacy/cbl/easy.html>.

If you want to create your own programs for LabPro, we encourage you to consult the Technical Reference document from the Vernier web site <http://www.vernier.com/mbl/labpro.html> for detailed explanations and additional information on the commands.


Sending Graphs and Data to Your Computer

Now that you have collected data, you can, of course, analyze data with your calculator. However, if you wish to analyze the data on a computer or print the data, you can connect the calculator to a computer and transfer the data from the calculator to the computer. To do this you will need software on the computer and a TI-GRAPH LINK cable.

You have three options when sending data back to your computer. One option is to simply use the TI-GRAPH LINK software for printing graphs and data lists. The TI InterActive!™ program includes a word processor with integrated math system, TI graphing calculator functionality, symbolic computer algebra system, web browser, data editor with spreadsheet, and calculator connectivity. Alternatively, you can use the Vernier Graphical Analysis program along with the TI-GRAPH LINK cable. Graphical Analysis software (Macintosh or Windows version) has a file option for importing data lists from all of the LabPro-compatible TI graphing calculators. The Graphical Analysis program is not required for data collection with LabPro. However, most people have found it is a great tool for directly moving data from the graphing calculator to the computer for graphing and printing. It is also very good for performing mathematical modeling and finding the relationships between variables. Also, as shown below, it produces high quality graphs.

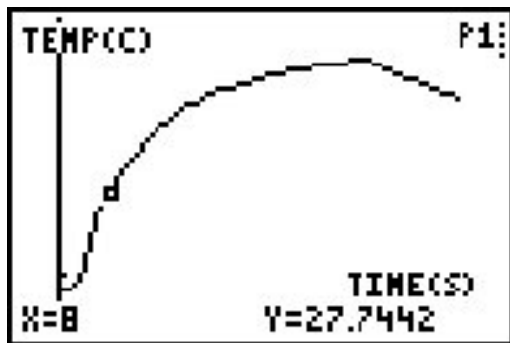

Using TI-GRAPH LINK software

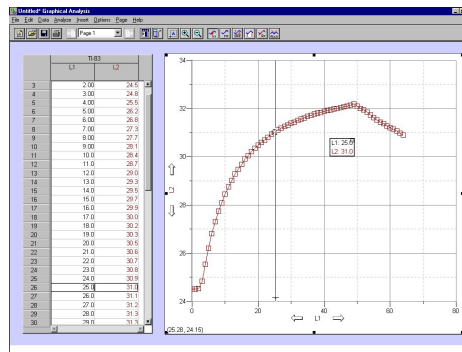



Using Graphical Analysis software


USING LOGGER PRO AND GRAPHICAL ANALYSIS SOFTWARE

You can use the Logger Pro or Graphical Analysis program to transfer the data on the calculator to your computer for analysis or printing. To do this, you need a TI-GRAPH LINK cable, which is part of the TI-GRAPH LINK package and is sold by Vernier Software & Technology and other Texas Instruments dealers. This cable connects the TI graphing calculator to the serial or USB port of your computer. Graphical Analysis 3.0 (order code GA). The purchase of this program includes a school or college department site license as well as a student license.

Note: Instructions for importing calculator data into Graphical Analysis is identical for Windows and Macintosh computers.

Importing procedures for TI-73, TI-73 Explorer, TI-83, TI-83 Plus, TI-83 Plus SE, TI-84 Plus, TI-84 Plus SE, TI-86, TI-89, TI-89 Titanium, TI-92, TI-92 Plus, Voyage 200:





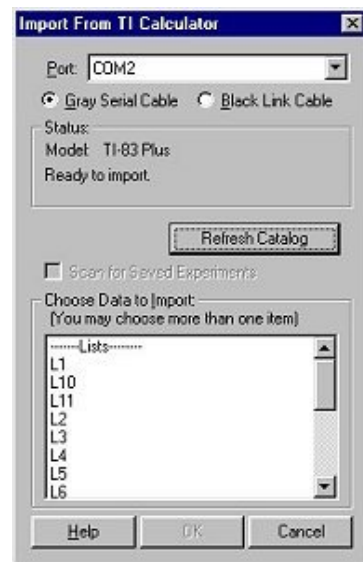

1. Connect the TI-GRAPH LINK cable to a free serial or USB port of the computer and to the TI calculator.
2. Turn on the calculator.
3. With Graphical Analysis running, pull down the File menu and choose Import From → TI Calculator or LabPro/CBL... or click the  button from the toolbar. The dialog box to the right should appear.
4. From the Port menu, choose the USB port or serial port (COM 1-4 on PC, modem or printer on Mac) to which the TI-GRAPH LINK cable is connected.
5. If you are using a PC serial cable, identify it is a gray or black cable.
6. Click on the Scan for Calculator button. The model you are using should now be identified, and you should see a message, “Ready to Import.”
7. Click on the data lists that you want to import. (To select more than one list on a Macintosh, hold down the  key while you click).
8. Click OK to send the data lists to the computer. The data will appear in columns in the data table. They will be labeled with the simple list names from the calculator. If you want to re-name them or add units, double-click on the column heading in the data table and enter new labels and units.




Importing procedures for the TI-82 and TI-85:

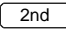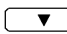

1. Perform steps 1-5 above.
2. Click on the Scan for Calculator button. The calculator model you are using should now be identified, and you should see a message “Waiting for you to transmit data from your calculator...”.
3. You can then choose which lists to transmit to the computer. Here’s how:

TI-82 Calculators:

On the TI calculator, press  [LINK], then select 2:Select All-... from the SEND menu. Use  to locate the lists on the SELECT menu. Position the arrow in front of a list you want to send to Graphical Analysis and press  to select it. More than one list may be selected in this manner. A  will appear beside each list to be sent. To deselect, press . The  will disappear.

Press , then select 1:Transmit. The lists will appear in the import window on the computer. They will be labeled with simple list names from the calculator.

TI-85 Calculators:

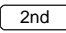On the TI calculator, press  [LINK], then press < SEND >. Press , then press <LIST>.

Use to locate a list you want to send and press < SELECT > to select it. More than one list may be selected in this manner. A ■ will appear beside each list that will be sent. To deselect, press . The ■ will disappear.

Press < XMIT > to transmit the list(s) to the computer. The lists will appear in import window. They will be labeled with simple list names from the calculator.

4. Click on the data list(s) that you want to import. (To select more than one list on a Macintosh, hold down the Apple key while you click).
5. Click OK to send the data lists to the computer. The data will appear in columns in the data table. They will be labeled with the simple list names from the calculator.

Appendix A: Battery and Adapter Information

OPERATING POWER REQUIREMENTS

- The LabPro is designed to operate with four AA alkaline batteries.
- Factors that affect battery life are the actual time that the LabPro is collecting data and the amount of current used by connected probes during your experiments.
- To extend battery life in the classroom we recommend that you use an approved power adapter.
- For long-term experiments out of the classroom, you can connect an external 6-Volt lantern battery to the LabPro. (See Connecting an External 6-Volt Battery on Page 40.)

WHEN TO REPLACE BATTERIES

The four AA batteries should be replaced when the low-battery message is displayed on the Main screen of DataMate. In addition, you can check the batteries at any time by choosing option CHECK BATTERY on the DataMate Tools screen.

Note: Save all collected data before removing batteries. All LabPro setup information and all collected data are lost if the batteries are removed and an optional AC adapter is not connected to the LabPro.

RECOMMENDED BATTERIES

- Four 1.5-Volt size AA alkaline batteries. Rechargeable batteries are not recommended.
- One 6-Volt lantern type. Recommended for long-term experiments performed outside the classroom that draw large amounts of current (for example, when using the Carbon Dioxide sensor). See Connecting an External 6-Volt Battery on Page 40 for instructions.

INSTALLING THE AA BATTERIES

- Holding the LabPro upright, push the latch on the battery cover down with your finger and pull the cover out.
- Replace all four AA alkaline batteries. Be sure to position them according to the diagram inside the battery compartment.
- Replace the cover.

CONNECTING AN OPTIONAL AC ADAPTER

- Connect one end of an approved adapter to the external power input connection located on the bottom left side of the LabPro.
- Plug the other end of the adapter into an electrical wall outlet.

APPROVED AC POWER ADAPTERS

The LabPro is designed to accept voltage input from an external AC-to-DC power adapter that provides a regulated 6 Volts DC output when plugged into an electrical wall outlet.

The model IPS power supply is an AC-to-DC power adapter approved for use with the LabPro. The TI-Presenter power supply can also be used with LabPro. The use of other power adapters is strongly discouraged.

To order an adapter, call Vernier at 1-503-888.837.6437 or visit the Vernier Web Store at <http://store.vernier.com>.

BUILDING AN EXTERNAL BATTERY ADAPTER CABLE

To build an external battery adapter cable, you will need a connector, 16-gauge wire (about 6 feet), and two alligator clips.

Note: At the time of this printing, the Radio Shack Coaxial DC Power Plug #274-1569 (5.5mm O.D., 2.1mm I.D.) or equivalent is an acceptable connector.

- Identify one 3-foot length of the wire as black (ground) and solder it to the insulated pin of the connector.
- Identify the other 3-foot length of wire as red and solder it to the outside of the connector.
- Connect an alligator clip to the open end of each wire.

CONNECTING AN EXTERNAL 6-VOLT BATTERY

- Connect one end of the external battery adapter to the external power input connection located on the right bottom edge of the LabPro.
- Connect the red lead to the positive (+) terminal of the battery. Connect the black lead to the negative (-) terminal of the battery.

Appendix B: DataMate and Calculator Specifics

There are functional and operational differences between the various types of TI graphing calculators. This Appendix describes those differences.

CALCULATOR DIFFERENCES

Because of the differences in memory between the calculators, there are some differences in functionality between the different versions of DataMate. See also the troubleshooting guide in Appendix E for information on the maximum number of data points each calculator can collect.

TI-73, TI-73 Explorer, TI-83 Plus, TI-83 Plus Silver Edition, TI-84 Plus, TI-84 Plus Silver Edition, TI-86, TI-89, TI-89 Titanium, TI-92, TI-92 Plus, and Voyage 200 versions support all of the DataMate functions.

TI-82

- The TI-82 version of DataMate supports only the following Auto-ID sensors: temperature, light, voltage, and CBR™ or the Vernier Motion Detector with pivoting head. Other sensors are supported, but due to memory constraints they are labeled as “Custom.” When selecting a custom sensor, you are responsible for entering the sensor’s slope and intercept. These values are listed with the booklet that’s included with the sensor, or can be found on the Vernier web site: <http://www.vernier.com/tech/manuals.html>.
- It supports all functions except SAVE/LOAD, SELECT REGION, and ADD MODEL.
- Supports up to 3 analog sensors.
- Supports the Motion Detector and only one analog sensor simultaneously.
- Adding auto-ID sensors increases the length of experiment and reduces the number of samples to save memory and keep experiment length the same.

TI-83

- The TI-83 version of DataMate supports all of the functions except SAVE/LOAD.
- DataMate will reduce the number of samples as needed to meet memory requirements.

TI-92

- DataMate will reduce the number of samples as needed to meet memory requirements.

Appendix C: LabPro Products, Calculators, and Sensors

Here are some of Vernier sensors that are available. Take a look at the Vernier catalog or check out the web site www.vernier.com for a full listing of LabPro and CBL 2 products.

SENSORS FOR THE LABPRO AND CBL 2

Sensor/Probe	Order Code
3-Axis Accelerometer	3D-BTA
25-g Accelerometer	ACC-BTA
Low-g Accelerometer	LGA-BTA
Barometer	BAR-BTA
Biology Gas Pressure Sensor	(use GPS-BTA)
CO ₂ Gas Sensor	C02-BTA
Colorimeter	COL-BTA
Conductivity Probe	CON-BTA
Current Probe System	DCP-BTA
Differential Voltage Probe	DVP-BTA
Digital Control Unit (<i>for LabPro & CBL 2</i>)	DCU-BTD
Digital Control Unit (<i>for original CBL</i>)	DCU-CBL (<i>CBL</i>)
Dissolved Oxygen Probe	DO-BTA
Drop Counter	VDC-BTD
Dual-Range Force Sensor	DFS-BTA
EKG Sensor	EKG-BTA
Exercise Heart Rate Monitor	EHR-BTA
Extra Long Temperature Probe	TPL-BTA
Gas Pressure Sensor	GPS-BTA
Heart Rate Monitor	HRM-BTA
Flow Rate Sensor	FLO-BTA
Instrumentation Amplifier	INA-BTA
Ammonium Ion-Selective Electrode	NH4-BTA
Calcium Ion-Selective Electrode	CA-BTA
Chloride Ion-Selective Electrode	CL-BTA
Nitrate Ion-Selective Electrode	NO3-BTA
Light Sensor	LS-BTA
Magnetic Field Sensor	MG-BTA
Microphone	MCA-BTA
Motion Detector	MD-BTD
O ₂ Gas Sensor	O2-BTA
ORP Sensor	ORP-BTA
pH Sensor	PH-BTA
Radiation Monitor	RM-BTD
Relative Humidity Sensor	RH-BTA
Respiration Monitor Belt (requires GPS-BTA or BGP-DIN)	RMB
Rotary Motion Sensor	RMS-BTD
Sound Level Meter	SLM-BTA
Stainless Steel Temperature Probe	TMP-BTA

Student Force Sensor	SFS-BTD
Student Radiation Monitor	SRM-BTD
Turbidity Sensor	TRB-BTA
Thermocouple	TCA-BTA
UVA Sensor	UVA-BTA
UVB Sensor	UVB-BTA
Vernier Photogate (Assembled)	VPG-BTD
Voltage Probe	VP-BTA

DO I NEED TO PURCHASE AN ADAPTER?

You do not need an adapter if you are purchasing new sensors for a LabPro or CBL 2. Simply buy the sensors with order codes that end in “-BTA” or “-BTD” and you’ll be ready to collect data. If you purchased older sensors, you may need an adapter. This chart should help. If you need additional help, call Vernier (888) 837-6437.

Sensor	Sensor Order Code	LabPro/CBL 2 Adapter	Original CBL Adapter
Analog (5-pin)	(---DIN)	DIN-BTA	DIN-BTA
Motion Detector (new, removable cable)	MD-ULI	MDC-BTD (Cable)	MDC-BTD (Cable),
Motion Detector (old, non-removable cable)	MD-ULI	CBL-RJ11	CBL-RJ11
Vernier Photogate	VPG-DG	PG-BTD (Cable)	CBL-P or CBL-2P
PASCO Photogate (has LED)	VPG-DG	DG-BTD	CBL-P or CBL-2P
PASCO older Photogate (no LED)	VPG-DG	Does Not Work	Does Not Work
Home-Made Photogate	PG-ULI	HMP-BTD	CBL-P or CBL-2P
Smart Pulley	ME97387	DG-BTD	CBL-P or CBL-2P
Radiation Monitor	RM-DG	DG-BTD	CBL-P
Student Radiation Monitor	SRM-DG	RAD-BTD (Cable)	CBL-P
Rotary Motion Sensor	RMS-ULI	ROT-BTD	Does Not Work

Appendix D: Support and Service Information

PRODUCT SUPPORT

Customers in the U.S., Canada, Puerto Rico, and the Virgin Islands

For general questions or technical questions contact Vernier.

phone: (503) 277-2299

e-mail: info@vernier.com

INTERNATIONAL ORDERS

All Vernier Software orders outside the US and Canada (except APO addresses) are handled by our international distributor, Ed. USA.

(Note that some Texas Instruments Dealers in other countries sell our products.)

Ed. USA, Inc.
5136 Calle Minorga
Siesta Key
Sarasota, FL 34242 USA
Fax: 941-349-2766
Phone: 941-349-1000
E-Mail: info@ed-usa.com

PRODUCT SERVICE

Contact Vernier before returning a product for service.

OTHER SERVICES

Visit the Vernier home page on the World Wide Web.

www.vernier.com

Visit the TI calculator home page on the World Wide Web.



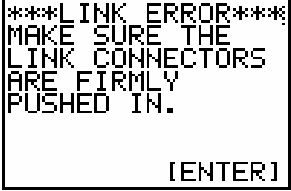
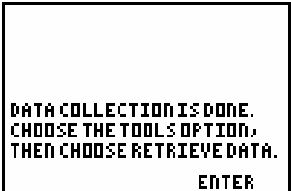
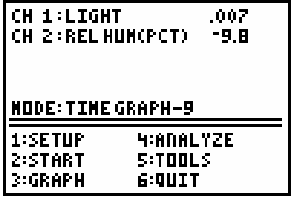
<http://education.ti.com/>





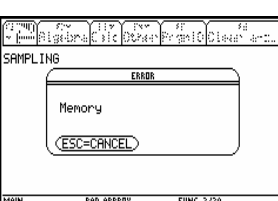
WARRANTY INFORMATION






Customers in the U.S. and Canada Only



All Vernier and Texas Instruments products carry a one-year limited warranty against defects in materials and workmanship. **In our 20 years, we have rarely charged a customer for repairs, except in cases of misuse.** All Texas Instruments products are covered by a one-year limited warranty against defects in materials or workmanship.

Appendix E: DataMate Troubleshooting Guide

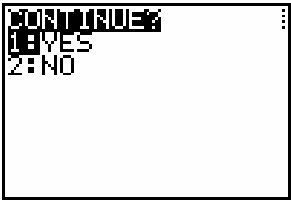Screen	Explanation
	<p>This screen appears when too much time has passed on a screen without activity. This timeout feature takes advantage of the Automatic Power Down (APD) feature of the calculator and LabPro to conserve battery power.</p> <ul style="list-style-type: none"> • Press YES to continue the program. • Press NO to quit.
 	<p>This screen appears when LabPro is not connected to the calculator or when LabPro needs new batteries.</p> <ul style="list-style-type: none"> • Check the connection between LabPro and your TI calculator. Firmly push in the link cable, and then choose 1: INTERFACE. • Check batteries in the LabPro. Disconnect the calculator from the LabPro; then press TRANSFER on the LabPro. If the LabPro does not make a sound or light the red LED, change the batteries in the LabPro. <p>If you choose 1: INTERFACE without correcting the problem, the link error screen appears.</p> <p>Check the connection and batteries as explained above and then press ENTER.</p> <p>If you choose 2: NO INTERFACE, then you can continue working in DataMate (viewing or analyzing previously collected data, etc.) but you won't be able to set up experiments or collect data.</p>
	<p>This screen appears when:</p> <ul style="list-style-type: none"> • The LabPro has collected data and that data has not been retrieved from the calculator. • Or • The user exits DataMate in the middle of data collection (possibly by pressing ON) and then restarts DataMate. <p>Press ENTER. Then choose one of the following:</p> <ul style="list-style-type: none"> • To retrieve the data, select TOOLS and select RETRIEVE DATA. • To delete the data, press <input type="button" value="CLEAR"/> to reset the LabPro.
	<p>The DataMate main screen shows a non-autoID sensor from a previous experiment even though the sensor is no longer connected. (For example, the screen on the left shows a Relative Humidity sensor even though the sensor was removed and DataMate was restarted.)</p> <p>Press <input type="button" value="CLEAR"/> to reset the LabPro to initial conditions. (In general, whenever you see something on the screen that does not look correct, press <input type="button" value="CLEAR"/> to reset.)</p>

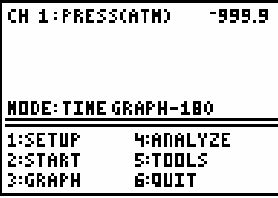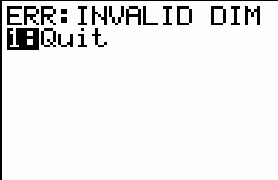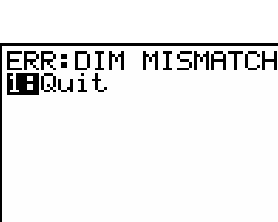Screen	Explanation																																				
	<p>This screen appears when the LabPro is disconnected from the calculator or when the LabPro loses power. When the LabPro and calculator are reconnected, the calculator may not recheck the sensor setup, and this error results.</p> <p>Press CLEAR to reset the LabPro to initial conditions (In general, whenever you see something on the screen that does not look correct, press CLEAR to reset.), and then set up the channel again. Or press GRAPH (On TI-89, TI-89 Titanium, TI-92, TI-92 Plus) to recalibrate channels but retain data collection parameters.</p>																																				
	<p>These four screens usually appear when there is not enough memory available in the calculator to collect any or all of the data and then graph it. Reduce the number of data points you are trying to collect.</p> <p>Following are <i>estimates</i> of the number of data points that can be collected if the calculator RAM memory has been reset prior to sending DataMate to the calculator:</p>																																				
	<table border="1"> <thead> <tr> <th>Calculator</th> <th>1 analog sensor</th> <th>2 analog sensors</th> <th>Digital sensor</th> </tr> </thead> <tbody> <tr> <td>TI-73</td> <td>~120</td> <td>~90</td> <td>~70</td> </tr> <tr> <td>TI-82</td> <td>98*</td> <td>98*</td> <td>98*</td> </tr> <tr> <td>TI-83</td> <td>~200</td> <td>~150</td> <td>~120</td> </tr> <tr> <td>TI-83 Plus</td> <td>998*</td> <td>~600</td> <td>~400</td> </tr> <tr> <td>TI-86</td> <td>~3000</td> <td>~2000</td> <td>~1500</td> </tr> <tr> <td>TI-89**</td> <td>~600/998***</td> <td>~600/998***</td> <td>~450/~800</td> </tr> <tr> <td>TI-92</td> <td>~300</td> <td>~200</td> <td>~150</td> </tr> <tr> <td>TI-92 Plus**</td> <td>~600/998***</td> <td>~600/998***</td> <td>~450/~800</td> </tr> </tbody> </table>	Calculator	1 analog sensor	2 analog sensors	Digital sensor	TI-73	~120	~90	~70	TI-82	98*	98*	98*	TI-83	~200	~150	~120	TI-83 Plus	998*	~600	~400	TI-86	~3000	~2000	~1500	TI-89**	~600/998***	~600/998***	~450/~800	TI-92	~300	~200	~150	TI-92 Plus**	~600/998***	~600/998***	~450/~800
Calculator	1 analog sensor	2 analog sensors	Digital sensor																																		
TI-73	~120	~90	~70																																		
TI-82	98*	98*	98*																																		
TI-83	~200	~150	~120																																		
TI-83 Plus	998*	~600	~400																																		
TI-86	~3000	~2000	~1500																																		
TI-89**	~600/998***	~600/998***	~450/~800																																		
TI-92	~300	~200	~150																																		
TI-92 Plus**	~600/998***	~600/998***	~450/~800																																		
	<p>* This is the limit of the TI-82/83 Plus calculator list.</p>																																				
	<p>** For the TI-89 and 92 Plus, the first value is the number of points that can be collected with 1.xx Calculator Operating System; the second value is the number that can be collected with 2.03 calculator operating system. The 2.03 Operating System optimizes the calculator memory; we recommend you upgrade to the 2.03 (or higher) code. The latest code is available at: http://www.ti.com/calc/docs/downloads.htm.</p> <p>*** This is the limit of the TI-89/92 Plus calculator data variable.</p> <p><i>Note: If you use DataMate's curve fit option, the memory available for collecting data will be less than that shown above because more calculator memory is required for curve fitting.</i></p> <p>Using a TI-89, TI-92 or TI-92 Plus, if the memory error occurs because too many data points were attempted, you must go into the memory management of the calculator and delete the "cbldata" data var. Then restart DataMate and begin data collection. Remember to reduce the number of points collected.</p>																																				

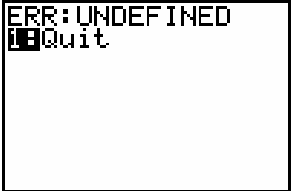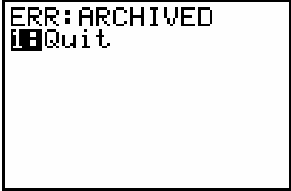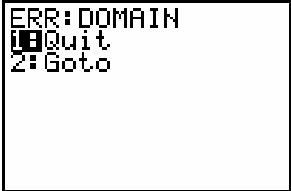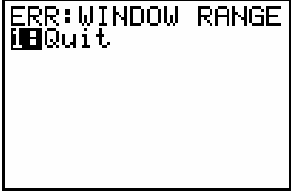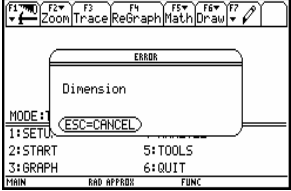Screen	Explanation
	<p>This screen usually appears when the user is running DataMate and one of the DataMate subprograms has been deleted from the calculator memory. (This applies to the TI-73, 82, 83, 86, 89, 93, and 92 Plus). All of the subprograms must be present in order for DataMate to function properly. (All related programs begin with "DATxxxx.")</p> <p>Reset the RAM on the calculator, then transfer the DataMate program from LabPro to the calculator and begin again.</p>
	<p>This screen appears on a TI-83 Plus calculator when one of the variables accessed by the DataMate app has been archived in the calculator memory. These variables should not be archived:</p> <ul style="list-style-type: none"> lists: L1 - L11, list C, list M real: A - Z matrix: [A] string: Str0 - Str6 <p>Go into Memory Management and unarchive any of the above variables.</p>
	<p>You have attempted a calculation outside the valid range. The most common cause of this error is trying to perform a power curve fit on Time Graph data. In Time Graph, DataMate collects a data point at time $x=0$. When the curve fit equation tries to divide by the 0, this error occurs.</p> <p>The easiest way to correct this is to use the SELECT REGION option to eliminate the $x=0$ point from the graph. Then try the power curve fit option again.</p>
	<p>The calculator tried to draw a graph but was not able to use the window settings. This problem can occur if you collect data and the data does not change (for example, the temperature does not change). If DataMate tries to autoscale the graph of this data (which it usually does), the calculator may not be able to set the y-axis scale.</p> <p>Select quit, then recollect the data. Make certain that the data changes for DataMate to graph correctly.</p>
	<p>This screen can appear when running DataMate on a TI-89/89 Titanium/92/92 Plus. It is caused by a loss of communications between the calculator and the LabPro and usually means there is a problem with the link port on the calculator.</p> <p>Check that the cable is securely connected to the calculator and the LabPro, then restart the program.</p>

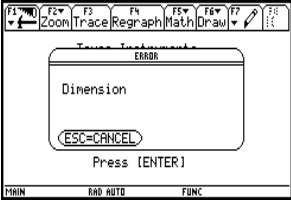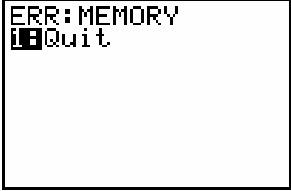Screen	Explanation
	<p>This screen appears when the user tries to run the Ranger program using the TI-89/89 Titanium/92/92 Plus after using the DataMate program.</p> <p>This is caused by a conflict in some information left over in List 5. The information in the list cannot be used properly, so the calculator gives a Dimension error. To correct this, go into the calculator's memory management and delete List 5 (L5).</p>
	<p>This screen can appear on the TI-83 Plus. It can be caused by running the DataMate app while the Interactive Graphing app is loaded and turned on.</p> <p>Turn off the Interactive Graphing app before running DataMate. Also, go into memory management and check the programs listing. There will be a program listed that has a strange character as its name. Reset the RAM on the calculator before doing anything else with the calculator.</p>
<p>When using the TI-82 with a motion detector and two other analog sensors, the data from the sensor in Channel 2 is not collected.</p>	<p>The TI-82 has the capacity for only six lists, so there are not enough lists available to collect data on all channels. When using a motion detector, you can only use one analog sensor in Channel 1.</p>
<p>The sensors and Time Graph Mode were set up in DataMate. Then Triggering was set up. When data collection started, the live graph did not display.</p>	<p>When triggering is selected, the LabPro does not allow a live graph.</p>

Appendix F: DataMate Sensor Setup Default Settings

The table that follows shows the default settings used by the DataMate program as well as the Quick Setup feature of LabPro. This sensor and calibration information applies only to Vernier and TI sensors. The use of other manufacturers' sensors may require calibration or the manual entry of that sensor's calibration information into DataMate.

Note: Default settings apply only when using a single sensor that is connected to Channel 1.

Sensor Name	Short Name	Y-Min	Y-Max	Sample Interval (in seconds)	No. Of Samples
Dir connect Temp (°C)	TEMP(C)	-15	110	1	180
Dir connect Temp (°F)	TEMP(F)	0	250	1	180
Extra Long Temp (°C)	TEMP(C)	-50	150	1	180
Stainless Temp (°C)	TEMP(C)	-20	125	1	180
Stainless Temp (°F)	TEMP(F)	-5	260	1	180
Thermocouple (°C)	TEMP(C)	-200	1400	1	180
pH	PH	0	14	2	60
Conduct 200 (µS)	CONDUCT (MICS)	0	200	1	180
Conduct 100(mg/l)	TDS(MG/L)	0	100	1	180
Conduct 2000(µS)	CONDUCT (MICS)	0	2000	1	180
Conduct 1000(mg/l)	TDS(MG/L)	0	1000	1	180
Conduct 20000(µS)	CONDUCT (MICS)	0	20000	1	180
Conduct 10000(mg/l)	TDS(MG/L)	0	10000	1	180
Gas Pressure (kPa)	PRESS (KPA)	50	150	10	90
Gas Pressure (mm Hg)	PRESS (MMHG)	400	1200	10	90
Gas Pressure (atm)	PRESS (ATM)	0.5	1.6	10	90
Gas Pressure (in Hg)	PRESS (INHG)	0	65	10	90
Pressure (kPa)	PRESS (KPA)	0	700	1	180

Sensor Name	Short Name	Y-Min	Y-Max	Sample Interval (in seconds)	No. Of Samples
Pressure (atm)	PRESS (ATM)	0	700	1	180
Pressure (mm Hg)	PRESS (MMHG)	0	5200	1	180
Bio Pressure (kPa)	PRESS (KPA)	76	156	10	90
Bio Pressure (mm Hg)	PRESS (MMHG)	550	1200	10	90
Bio Pressure (atm)	PRESS (ATM)	0.75	1.6	10	90
Bio Pressure (in Hg)	PRESS (INHG)	20	50	10	90
Dual R Force (5N)	FORCE(N)	-5	5	0.05	100
Dual R Force (10N)	FORCE(N)	-10	10	0.05	100
Dual R Force (50N)	FORCE(N)	-50	50	0.05	100
Student Force (N)	FORCE(N)	-40	10	0.05	100
EX Heart Rate (bpm)	Heart RT(BPM)	45	170	5	180
Heart Rate (bpm)	Heart RT(BPM)	45	170	5	180
25G Accel (m/s ²)	ACCEL (M/S ²)	-250	250	0.05	100
Low G Accel (m/s ²)	ACCEL (M/S ²)	-50	50	0.05	100
Colorimeter	ABSOR-BANCE	0	0.6	5	180
CO ₂ Gas (ppm)	CO ₂ GAS (PPM)	0	5000	10	30
CO ₂ Gas (ppt)	CO ₂ GAS (PPT)	0	5	10	30
CBL Microphone	MICRO-PHONE	0	5	1.00E-04	200
ULI Microphone	MICRO-PHONE	0	5	1.00E-04	200
MPLI Microphone	MICRO-PHONE	-5	5	1.00E-04	200
TI Light Sensor	LIGHT	0	1	0.05	180

Sensor Name	Short Name	Y-Min	Y-Max	Sample Interval (in seconds)	No. Of Samples
Light 600(lx)	LIGHT(LX)	0	600	0.05	180
Light 6000(lx)	LIGHT(LX)	0	6000	0.05	180
Light 150000(lx)	LIGHT(LX)	0	150000	0.05	180
D. Oxygen (mg/l)	DO(MG/L)	4	12	2	60
EKG	EKG	-0.5	4	0.01	200
CA ISE (mg/l)	CA(MG/L)	0	40000	1	180
NH4 ISE (mg/l)	NH4(MG/L)	0	18000	1	180
NO3 ISE (mg/l)	NO3(MG/L)	0	14000	1	180
CL ISE (mg/l)	CL(MG/L)	0	36000	1	180
Flow Rate (m/s)	FLOW RT (M/S)	0	4	1	180
Flow Rate (ft/s)	FLOW RT (FT/S)	0	13	1	180
Respiration (bpm)	RESP RT (BPM)	0	30	10	180
Turbidity (ntu)	TURBID (NTU)	0	50	1	180
C V Current (A)	CURRENT (A)	-0.6	0.6	0.1	180
C V Voltage (V)	VOLTAGE (V)	-6	6	0.1	180
Voltage -10 to +10 (V)	VOLTAGE (V)	-10	10	0.1	180
Voltage 0 to 5 (V)	VOLTAGE (V)	0	5	0.1	180
Hi Magnet Fld (mT)	MAGNET F(MT)	-0.32	0.32	0.05	180
Hi Magnet Fld (g)	MAGNET F(G)	-3.2	3.2	0.05	180
Lo Magnet Fld (mT)	MAGNET F(MT)	-10	5	0.05	180
Lo Magnet Fld (g)	MAGNET F(G)	-100	50	0.05	180
Barometer (kPa)	BARO(KPA)	80	110	600	180
Barometer (mm Hg)	BARO (MMHG)	600	800	600	180

Sensor Name	Short Name	Y-Min	Y-Max	Sample Interval (in seconds)	No. Of Samples
Barometer (in Hg)	BARO (INHG)	24	32	600	180
Barometer (mBar)	BARO (MBAR)	810	1060	600	180
Relative Humidity (%)	REL HUM (PCT)	0	100	600	180
Oxygen Gas (%)	O2 GAS (PCT)	15	25	15	40
Oxygen Gas (%)	O2 GAS (PPT)	150	250	15	40
Custom 0 to 5 (V)	CUSTOM	0	5	1	180
Custom -10 to +10 (V)	CUSTOM	-10	10	1	180
Motion (m)	MOTION (M)	0	6	0.05	100
Motion (ft)	MOTION (FT)	1	20	0.05	100
Current Probe (A)	CURRENT (A)	-10	10	0.1	180
Resistance (ohm)	RES (OHM)	0	100000	0.1	180
Sound Level	Sound(DB)	0	130	6	100

The values in the table are valid for all calculator versions of DataMate except the TI-82, TI-83, and TI-92. Because of memory limitations on these calculators, the number of data points is reduced.

- For the TI-82, all default number of samples will be limited to 99 points.
- For the TI-83, the number of samples will be reduced if it exceeds 200 points and multiple channels are set up.