

## LabQuest Viewer Release Notes

Version 2.0/2.0.1

March, 2015

### System Requirements

The LabQuest Viewer software is compatible with the following computer platforms:

- Version 2.0 – Macintosh OS X 10.8, 10.9, and 10.10
- Version 2.0.1 – Windows 7 SP1, Windows 8, and Windows 8.1

LabQuest Viewer 2.0/2.0.1 requires approximately 5.0/3.0 MB (Mac/Windows) disk storage; you need less than 10 MB free space to allow the installation to complete successfully.

Computer/projector display resolution 1024x768 or higher is recommended. The minimum display resolution is 800x600.

### Install Notes

The LabQuest Viewer 2.0/2.0.1 software makes use of Apple's Bonjour software for automatic device discovery. While Windows computer software can be used without Bonjour, requiring the user to manually identify devices for connection, it is not recommended.

### What's Changed in LabQuest Viewer 2.0.1 (*Windows only*)

- Fixed issue where LabQuest Viewer could leave LabQuest unresponsive to taps.

### What's Changed in LabQuest Viewer 2.0

#### ***New Features***

- New My Lab view displays live images of all LabQuests in your lab.
- My Lab Setup allows you to add, arrange, and remove LabQuest devices in My Lab.
- Persistence of My Lab setup and software size/position values between sessions.
- Previous/Next buttons support quickly moving through LabQuest units while presenting.
- Select up to four LabQuest for side-by-side screen comparison.

#### ***Other Changes***

- Full software support for Mac OS X 10.10.

Version 2.0 does not support Windows XP or Vista. Users of these operating systems must continue using LabQuest Viewer software version 1.1.